



# ODYSSEY<sup>®</sup>

## Tutorial Manual

Version 1.0



**LI-COR<sup>®</sup>**  
Biosciences

## **Notice**

The information contained in this document is subject to change without notice. LI-COR MAKES NO WARRANTY OF ANY KIND WITH REGARD TO THIS MATERIAL, INCLUDING, BUT NOT LIMITED TO THE IMPLIED WARRANTIES OF MERCHANTABILITY AND FITNESS FOR A PARTICULAR PURPOSE. LI-COR shall not be liable for errors contained herein or for incidental or consequential damages in connection with the furnishing, performance, or use of this material.

This document contains proprietary information that is protected by copyright. All rights are reserved. No part of this document may be photocopied, reproduced, or translated to another language without prior written consent of LI-COR, Inc.

## **Printing History**

Publication Number 984-11074

June, 2010

LI-COR, Odyssey, MPX and IRDye trademarks contained in the Software Product are trademarks or registered trademarks of LI-COR, Inc. Third party trademarks, trade names, and product names may be trademarks or registered trademarks of their respective owners. You may not remove or alter any trademark, trade names, product names, logo, copyright or other proprietary notices, legends, symbols, or labels in the Software Product. This EULA does not authorize you to use LI-COR's or its licensors' names or any of their respective trademarks.

LI-COR is an ISO9001 registered company. ©2010 LI-COR Inc. All rights reserved. Specifications subject to change. LI-COR, Odyssey, FieldBrite, MPX and IRDye are trademarks or registered trademarks of LI-COR, Inc. Odyssey Fc Imager, FieldBrite, and IRDye reagents are covered by U.S. patents, foreign equivalents, and patents pending.

# Table of Contents

---

## Table of Contents

<b>Chapter 1: Getting Started</b>	<b>1</b>
Installing the Software	1
Setting up a Work Area	1
Connecting to the Odyssey® Fc	1
Using Key Tips	2
<b>Chapter 2: System Overview</b>	<b>3</b>
Odyssey Fc System: Near-Infrared Fluorescence and Chemiluminescence	3
Odyssey Fc Detection System	3
Table Format	3
Starting Scans and Analyzing Images	4
What's Next?	4
<b>Chapter 3: Acquiring an Image</b>	<b>5</b>
Inserting the Sample	5
Acquiring the Image	5
Viewing the Image and Data	7
<b>Chapter 4: Importing an Image</b>	<b>9</b>
Importing an Image	9
<b>Chapter 5: Manipulating an Image</b>	<b>11</b>
Viewing the Image	11
Adjusting the Look of the Image	11
Rotating the Image	14
Cropping the Image	15
<b>Chapter 6: Analyzing a Blot or Gel</b>	<b>17</b>
Adding and Manipulating Shapes Manually	17
Using the Profiles	20
Subtracting the Background	20
Annotating the Image	21

<b>Chapter 7: Performing Western and MPX™ Western Analyses.....</b>	<b>23</b>
Applying a Western Analysis.....	23
Setting the Boundary .....	24
Setting the Lanes.....	24
Setting the Marker .....	26
Finding the Bands Automatically .....	27
Editing Bands Manually.....	28
Subtracting the Background.....	29
Reviewing the Lanes Table.....	29
Reviewing the Bands Table .....	30
Applying the MPX Western Analysis.....	30
Reviewing the MPX Lanes Table .....	32
<b>Chapter 8: Working with the Tables.....</b>	<b>33</b>
Working in the Images Table .....	33
Adding or Deleting a Column .....	34
Editing the Image Title .....	35
Adding Custom Columns.....	35
Pasting Data in a Column .....	36
Copying Data to a Spreadsheet.....	37
Sorting by Column.....	37
Filtering the Table Manually .....	38
Defining Filters .....	38
Viewing the Shapes Table .....	39
<b>Chapter 9: Exporting Data, Printing Images, and Generating Reports.....</b>	<b>41</b>
Exporting Acquisition Data and Images .....	41
Printing Images .....	42
Generating Reports.....	43

# Chapter 1: Getting Started

---


## Installing the Software

Install the Image Studio Software and the Image Studio Odyssey® Fc Acquisition Key first. For added features, install the Image Studio Western Analysis Key. Once the software is installed, double-click the shortcut that appears on the Desktop to start the Image Studio software.

## Setting up a Work Area


When starting the software for the first time, you will need to create the first Work Area. The Work Area is a folder on your hard drive or network where all of your images, analyses and settings will be stored. Click on **Browse...** in the *Set Active Work Area* menu to open a browse window where you can select a folder on your hard drive or network to use as your Work Area. This folder will now appear in the *Available Work Areas* window every time you open the software. Select the folder and click **OK** to set it as your Work Area. (To delete a Work Area, select the folder in the *Available Work Areas* window and press the *Delete* key on your keyboard.)

**Note:** Each user should create their own Work Area, as the settings for the instrument, analysis, image display, etc. from the last session will be applied to the next session.

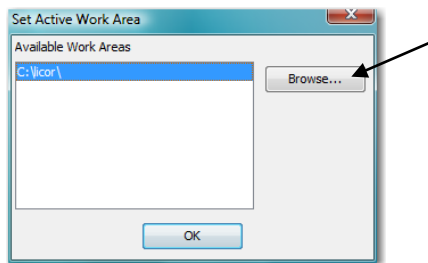


## Connecting to the Odyssey Fc

Once the software starts, it searches for and discovers the Odyssey Fc Imager on the same LAN automatically. If you experience connection problems, be sure that IPv6 is enabled on the computer. To manage the status of the instrument, click on the *Image Studio Application* button and hover over **Instrument**.

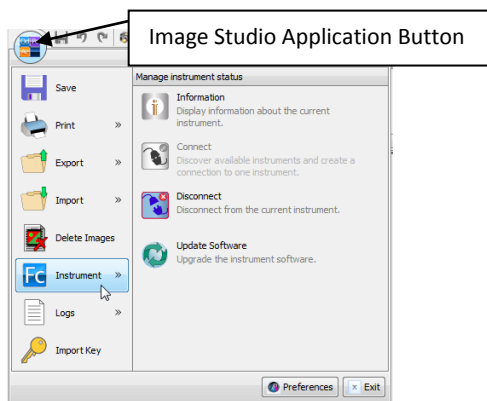


## Using Key Tips

The keyboard can be used to access the tabs and other features of the Image Studio software by activating Key Tips. Press **ALT** to display the key tip badges for all ribbon tabs and the application quick launch buttons. Press the letter key next to the application quick launch buttons to select that application.

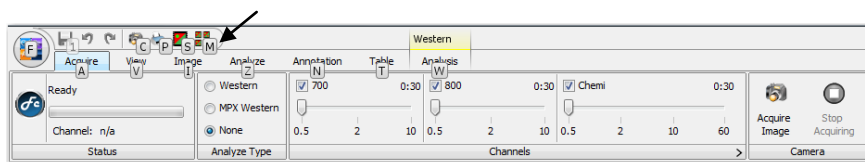

Press the letter key next to the ribbon tabs to view the ribbon for the selected tab and the key tips for the commands on that ribbon. Press the letter key next to the commands to activate the commands or open another menu.

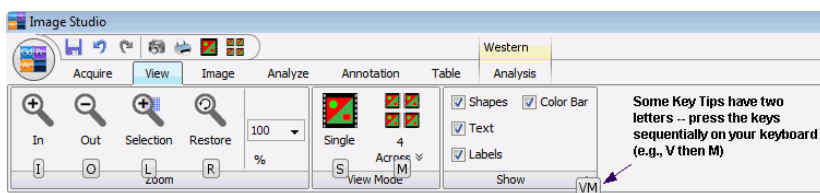

In this tutorial, directions will be given to select the tabs and commands by clicking with the mouse. Press **ALT** at any time to use the Key Tips to select the tabs and commands with the keyboard.

# Chapter 2: System Overview

---

## **Odyssey® Fc System: Near-Infrared Fluorescence and Chemiluminescence**

The Odyssey Fc System detects IRDye® Infrared Dyes at independent wavelengths for two-color detection and quantification in the infrared. Chemiluminescent substrates can also be detected and a qualitative image acquired. Each channel is acquired sequentially. An example of the use of the infrared channels with the chemiluminescent channel is imaging a marker lane stained with Coomassie in the 700 nm channel combined with the qualitative chemiluminescent image.

The latest LI-COR protocols are posted on-line at <http://biosupport.licor.com>. Files are in Adobe® Acrobat® format (\*.PDF), requiring a copy of Adobe Reader® to open the files. Adobe Reader is a free download at <http://www.adobe.com>.

## **Odyssey Fc Detection System**

The Odyssey Fc laser module contains a 685 nm and a 785 nm laser. Image acquisition by the LI-COR CCD camera consists of images for two fluorescent probes in the 700 and 800 nm channels, as well as an image for chemiluminescent substrates in the Chemi channel (assuming all channels are enabled). In the software, acquisition of the three channels is sequential and can be selected separately or in any combination.

The quality of Odyssey Fc images is enhanced by a patented filtering system that dramatically reduces noise before detection by the CCD detector. Infrared signal detection has also been optimized for IRDye Infrared Dyes, which eliminates the need for filter selection by the user before imaging. The unique proprietary FieldBrite™ XT technology used in the Odyssey Fc Imager and developed exclusively by LI-COR acquires images without saturated pixels on the first attempt. More than six logs (22 bits) of dynamic range are available for each image with no user adjustments.

## **Table Format**

Images acquired are referenced in a table with multiple columns that can be used to filter, sort, and select the desired images. There are a number of column descriptions which allow filtering by such things as Acquire Time, Project, or Channels. There are also three custom field columns. Images can be selected for analysis, printing, or export. Data from an analysis of the image are stored in tables with multiple descriptive columns as well. Image Studio has several options for exporting image and table data, and for generating reports.

## Starting Scans and Analyzing Images

For typical scans, a blot or gel is placed on the imaging tray and inserted in the imaging drawer. The scan is started from the computer using Image Studio software. In the software, the desired channel or channels are selected, along with the length of the integration time for each channel. At the end of a scan, image data are referenced in the Images Table with multiple descriptive columns.

When a scan is complete, you can open the *View* ribbon to modify how the image looks on the screen. Use the *Image* ribbon to create a modified copy of the image. In the *Analyze* ribbon, shapes are manually added and used for quantitation. The *Annotation* ribbon is used to add text and arrows to the image. Installing the Western Key adds the *Western Analysis* and *MPX Analysis* ribbons. These applications have enhanced features for adding lanes, bands, and markers to a Western or MPX™ Western blot.

## What's Next?

In Chapters 3 and 4, you will learn how to acquire and import images. Manipulating and analyzing an image is covered in Chapters 5 and 6. The Western Key application is discussed in Chapter 7, followed by instructions on working with the Tables in Chapter 8. In the final chapter, you will export and print the image and data, as well as generate a report.



# Chapter 3: Acquiring an Image

---

*Before starting, check that the instrument is turned on and that the Image Studio software with the Western Key is running and connected to the instrument.*

## Inserting the Sample

The first step is to insert a sample into the instrument. For the Tutorial, you will image the field installation target.

- 1) Open the imaging drawer by pressing the Imaging Drawer open/close button.
- 2) For this exercise, place the field installation target that comes with the Odyssey® Fc instrument in the imaging drawer by aligning the notches in the target with the guide pins in the drawer.


**Note:** Do not use the imaging tray for placing the field installation target in the instrument.

- 3) Press the Imaging Drawer open/close button again to close the drawer.


**Note:** Techniques for placing Western blots or gels in an imaging tray to insert in the instrument are given in the Odyssey Fc Operator's Manual.

## Acquiring the Image


- 1) To acquire an image, first click on the **Acquire** tab to show the *Acquire* ribbon.
- 2) Hover over the first group to learn the status of the instrument. Right now, it should say *Ready*, indicating the instrument is connected and not currently acquiring an image.

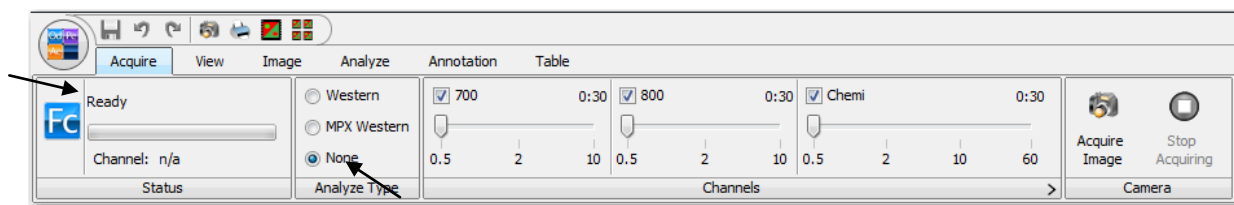


- 3) In the *Analyze Type* group, select **None** for this exercise.
- 4) In the *Channels* group, click the check boxes beside each channel (700, 800, and Chemi) so that a check mark appears, indicating each channel is selected for image acquisition.

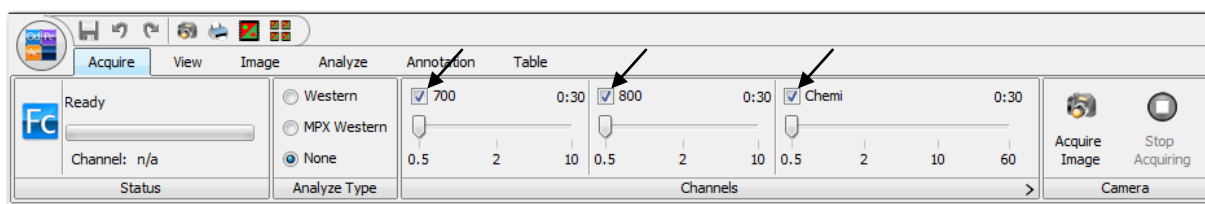

- 5) Select 0.5 minutes as the Integration Time for each channel by dragging the slider to 0.5 in each channel.

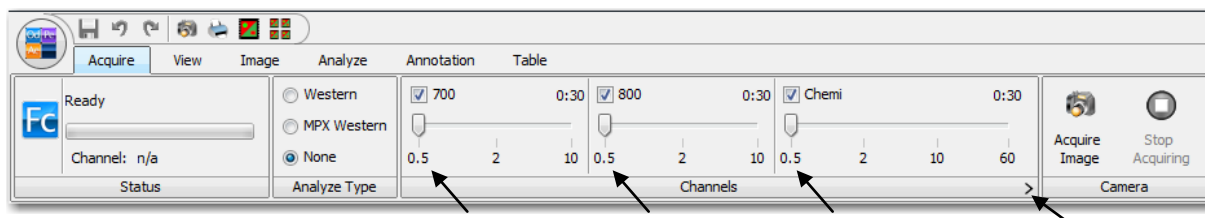


**Note:** In the Standard mode, only the shown times can be selected. To select an integration time that is not shown, click on the Expand Button at the end of the Channels bar to open the **Integration Time** menu. Click on **Extended** to change the Sliders Mode. Click and drag each Slider to select specific Integration Times.

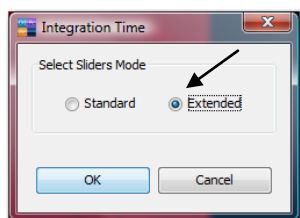

**Note:** Increasing the Integration Time will decrease the background noise and increase the signal-to-noise ratio.

- 6) Once the parameters have been set, click on **Acquire Image** to start the acquisition.

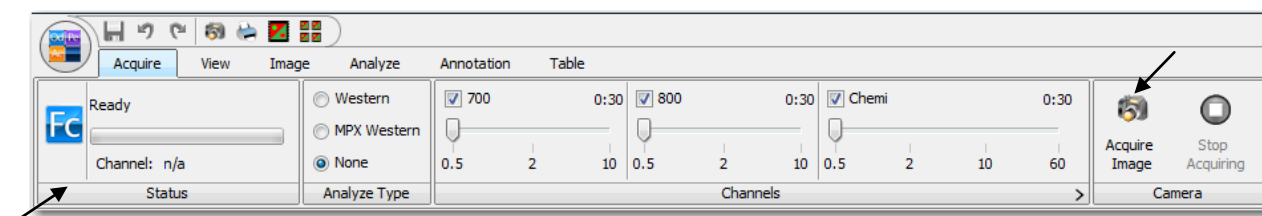


**Note:** The instrument acquires each channel separately. The Status group provides information on the imaging process.

- 7) If the acquisition needs to be stopped, click on **Stop Acquiring**. All existing and pending channel images are discarded.

## Viewing the Image and Data

When the acquisition is complete, the image will appear on the screen. The image data will appear in the table below in the Images Table. Change the Image Name to 'Tutorial' to easily find the image later.

- 1) Double-click on the cell under **Image Name** so that a cursor appears.
- 2) Key in 'Field Target' where the cursor appears. Press *Enter* on the keyboard when complete.
- 3) To view the image as it appears here,
  - a. Drag the middle dot in the Image LUTs to set K equal to 0 for the 700 and Chemi channels.
  - b. Adjust the max and min of the 700 channel by dragging the left and right dots to include all the pixels.
  - c. Refer to [Chapter 5: Manipulating an Image](#) for complete information on adjusting the look of the image.

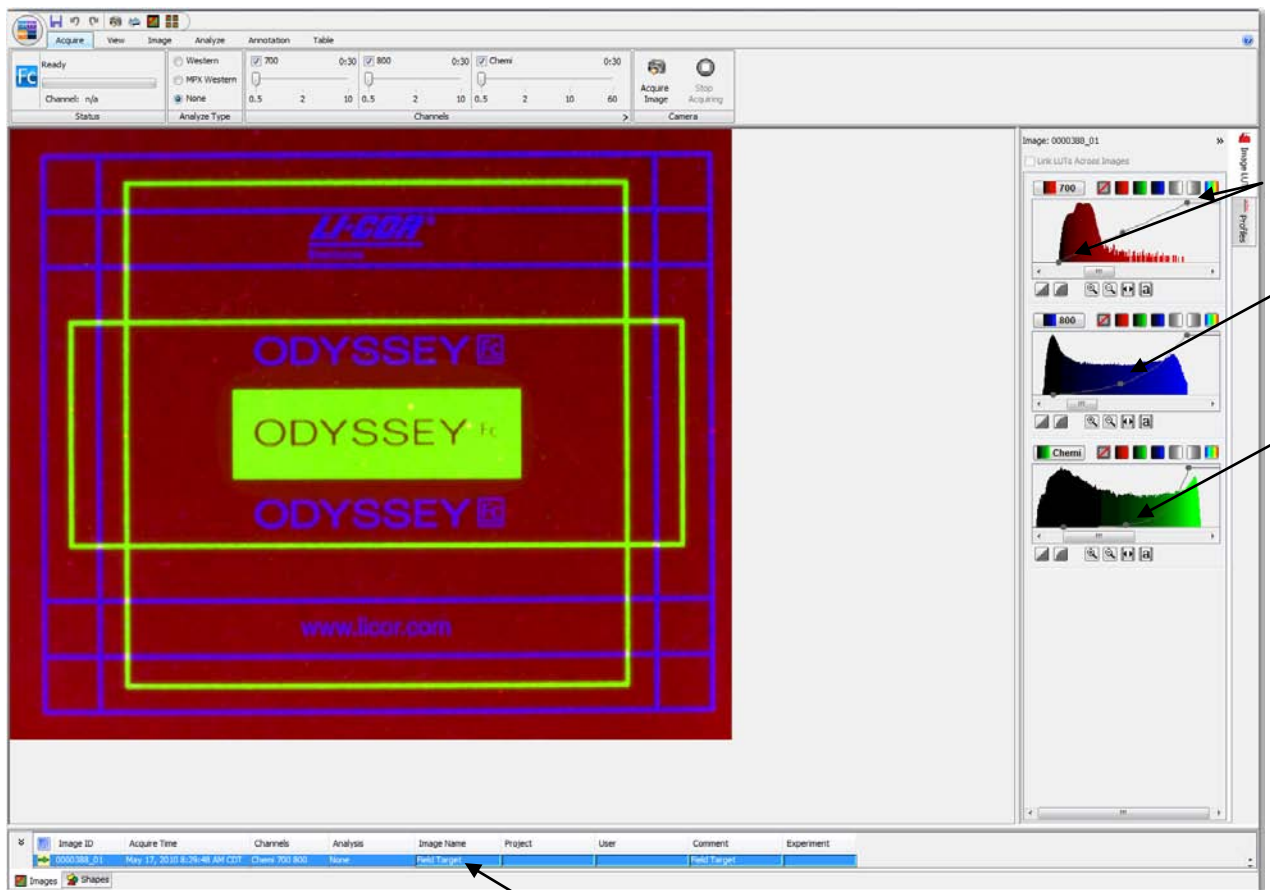



# Chapter 4: Importing an Image


---

*In Chapter 3, an image of the field installation target was acquired instead of a real Western blot or a gel for the purpose of the Tutorial. In this chapter, an image of a Western blot will be imported into your Work Area to use as an example for the next chapters.*

## Importing an Image


*The image needed for this tutorial is located on the Image Studio CD in a directory called Image Examples. The file name is '9999999\_01'.*

- 1) To import the image, click on the Application Menu Button and move the mouse to hover over **Import**. Click on **Copy From (Import Acquisition)** in the Import menu.

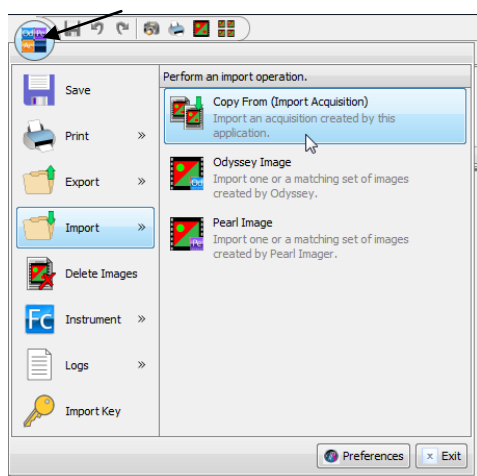


**Note:** Images from the Odyssey Imager can be imported into Image Studio software for analysis by clicking on **Import > Odyssey Image**.

- 2) Click once on the folder '9999999\_01' to select it. Click **Open** to display the image.

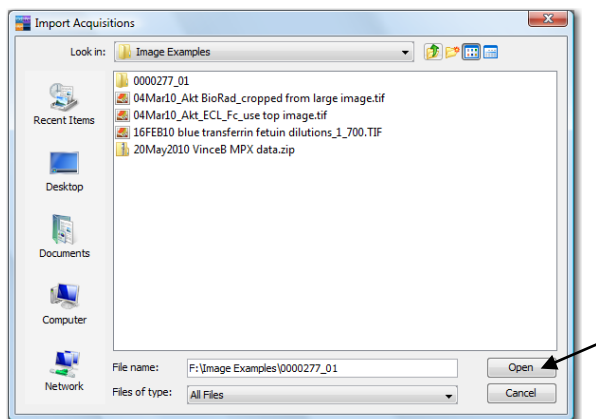


3) The image will open in the window and the file will appear in the Table.

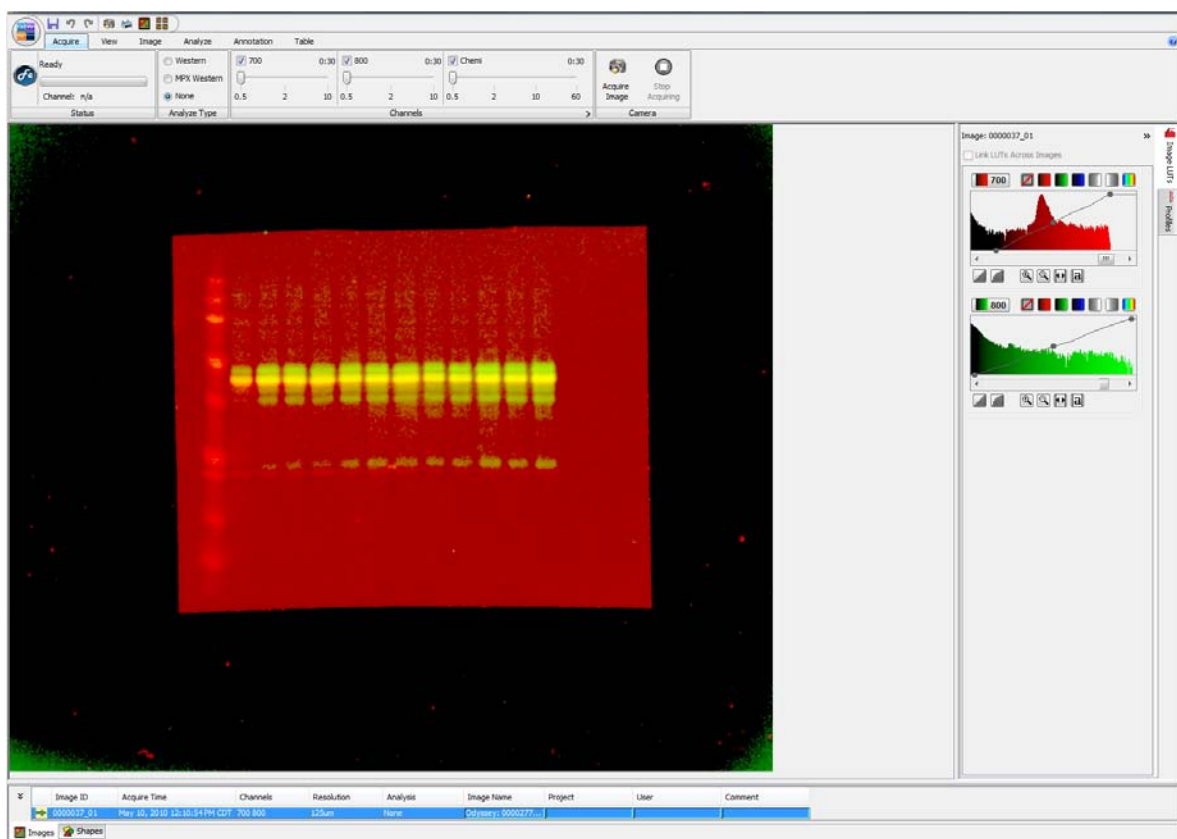

# Chapter 5: Manipulating an Image

## Viewing the Image


- 1) Click on the **View** ribbon tab to show the *View* ribbon.

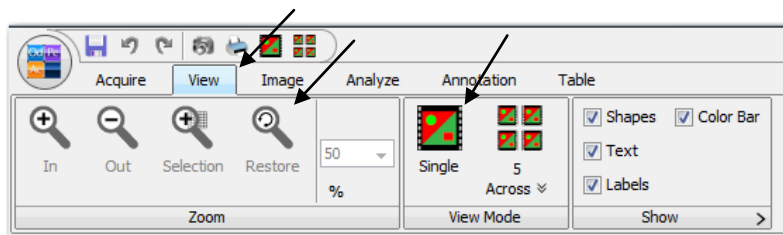

- 2) To increase and decrease the size of the image on the screen, click on the zoom icons or scroll the mouse wheel. For the tutorial, click on **Restore** to automatically fit the image to the window.
- 3) Select **Single** in the *View Mode* group to view and analyze a single image.

**Note:** Analysis cannot be done if the **Multiple Images** icon is selected.


- 4) In the *Show* group, click on all the boxes to show a check mark, indicating all items will appear on the image. To specify the location of the labels, click on the Expand button to open the *Image View Label Locations* menu.

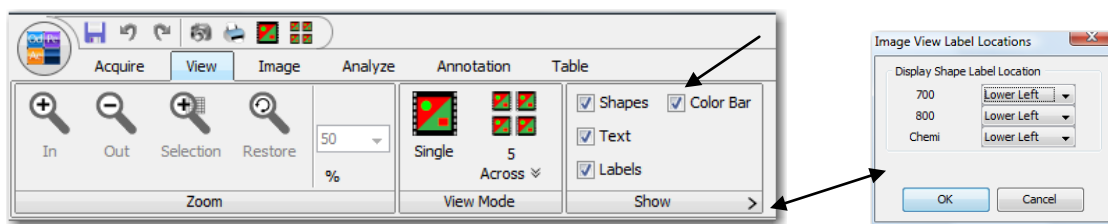

## Adjusting the Look of the Image


The *Look Up Tables* or *LUTs* show the histograms of the pixel intensities for each channel acquired for the image. The *LUTs* are on the right side of the screen.

- 1) Click once on the **Image LUTs** tab to view the histograms. To hide this view, click once on the Expand button.

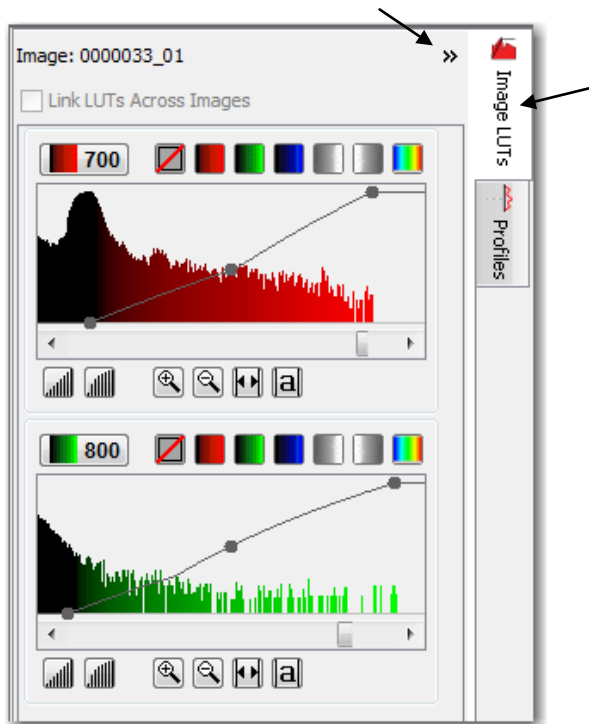

- 2) For the tutorial, the default red and green color scheme will be used. To select a different color scheme for each channel or to disable a specific channel, click once on the appropriate colored button above the graph.


**Note:** Multiple channels can be viewed, printed, and exported in gray scale.

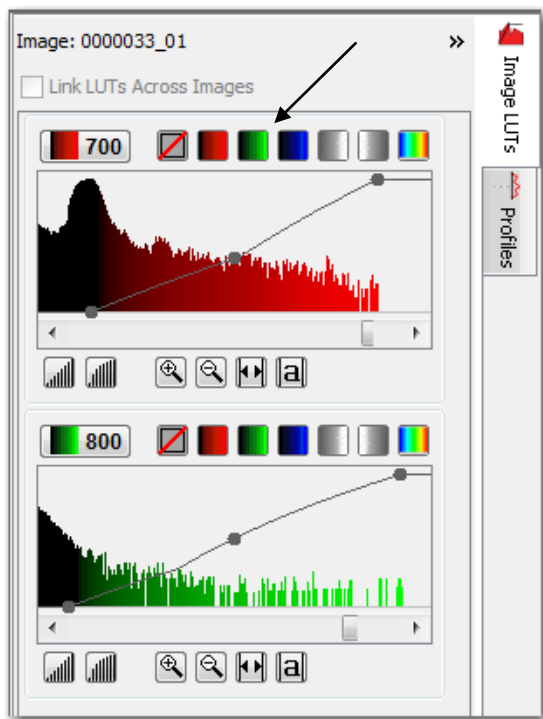



- 3) To adjust the intensity of the image,
  - a. Click the *Auto Adjust* button under each graph to automatically set the min and max points of the display mapping.
  - b. Click and drag the left (min) dot on each graph to adjust the lower threshold of pixel values that will shade to the same color on the viewed image. Shading the lower intensity pixels to the same color will visually create a cleaner background on the image.


**Note:** Changing the visual image will not change the data collected for that image or the quantification of that data. It only changes the mapping of raw image pixels to display pixels.

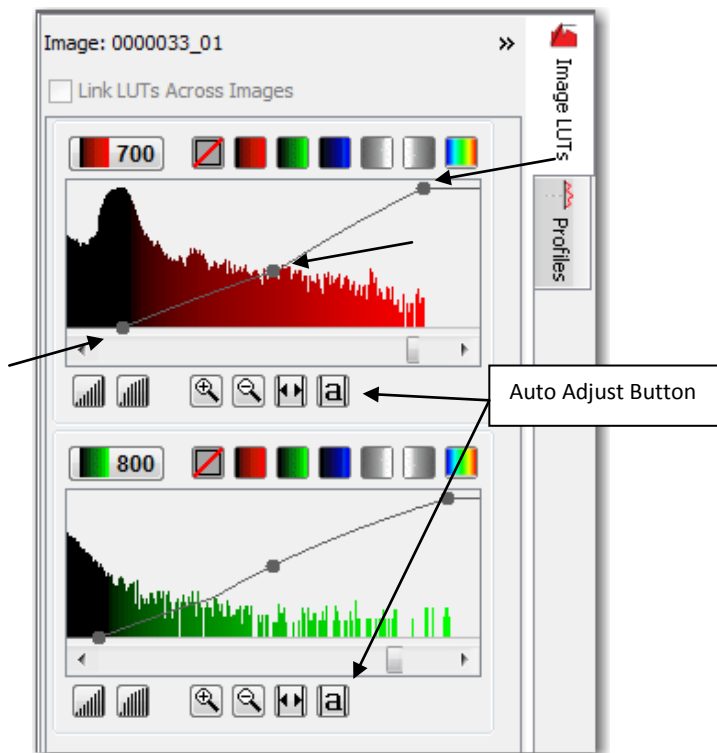

- c. The right (max) dot on each graph can be left at the auto adjusted value. It adjusts the upper threshold of pixel values that will shade to the same color on the viewed image.
  - d. Click and drag the middle dot on each graph vertically to change the K value.


**Note:** Further explanation of the K value can be found in the on-line Help system by clicking the question mark in the upper right hand corner of the screen.

The image should now look similar to this one.

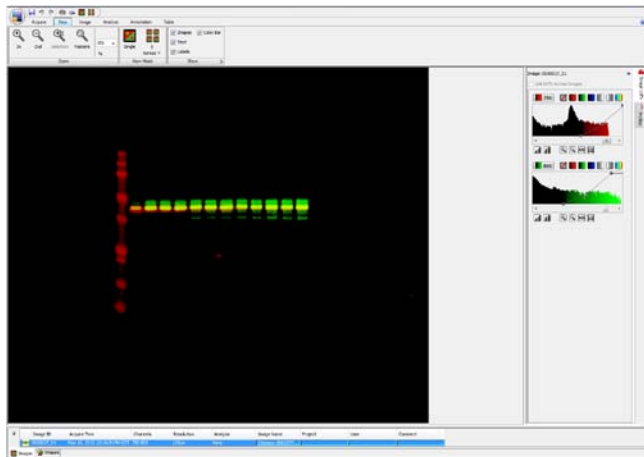


## Rotating the Image

- 1) Click on the **Image** ribbon tab to show the *Image* ribbon.

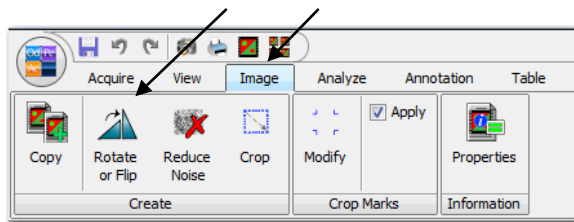

- 2) Click on the **Rotate or Flip** icon to display the *Rotate or Flip Options* menu.

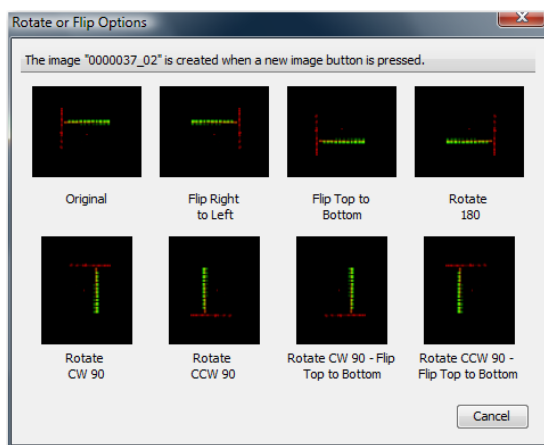


- 3) Select **Original** by moving the mouse to the thumbnail and clicking once on the highlighted box.

**Note:** A new image is created in the desired format and a new file for that image appears in the Table.


## Cropping the Image

- 1) To crop the image, first click on the box next to **Apply** so that a check mark appears in the check box. Small marks indicating the corners of the cropped area appear on the image.

**Note:** These small marks will not appear on the image when it is in gray scale. For cropping, change the color of the image.

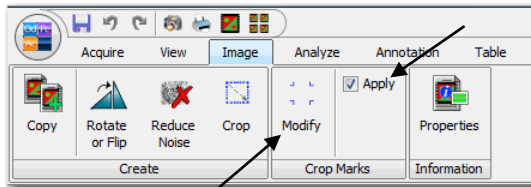


- 2) Next, click on **Modify** to open the *Edit Image Crop Marks* menu. Two boxes appear on the image in dotted lines.

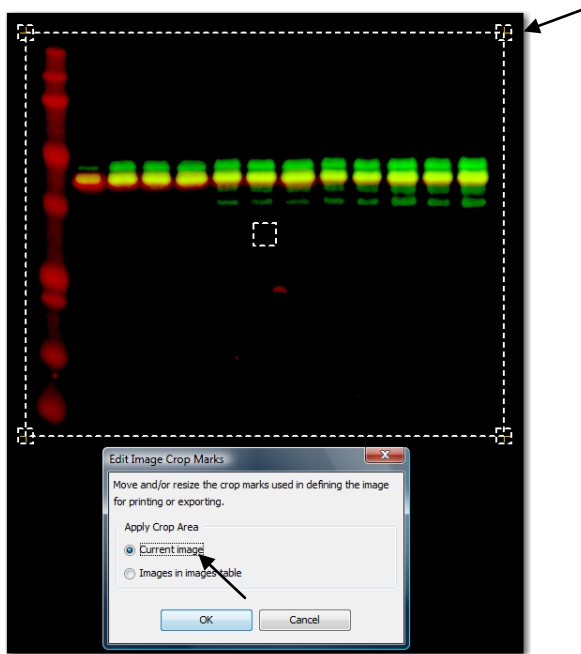


- 3) Click and drag the small box in the middle of the image to center the crop marks. Click and drag the corners of the large box to surround all of the marker and bands.
- 4) Click on **Current Image** to mark this image with these crop marks and click **OK**. Small crop marks indicating the corners of the cropped area appear on the image at the new location.

**Note:** If **Images in images table** is selected, this crop area is applied to all images in the *Images Table*.

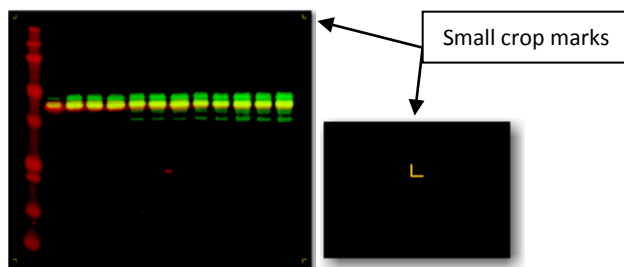


**Note:** These small crop marks will appear on the image as long as **Apply** has a check mark by it. The crop area will be applied to any print or export action. To remove the small crop marks and print or export the full image, click **Apply** to remove the check mark.

5) To create a new image of the cropped area, click **Crop**.

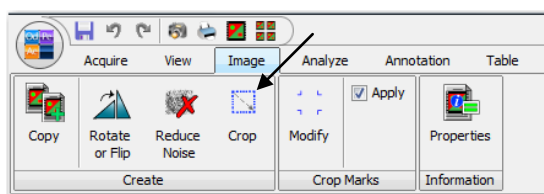


6) The cropped image appears in the window and a new file for that image appears in the Table. Click **OK** in the *Crop Image Completed Successfully* box.

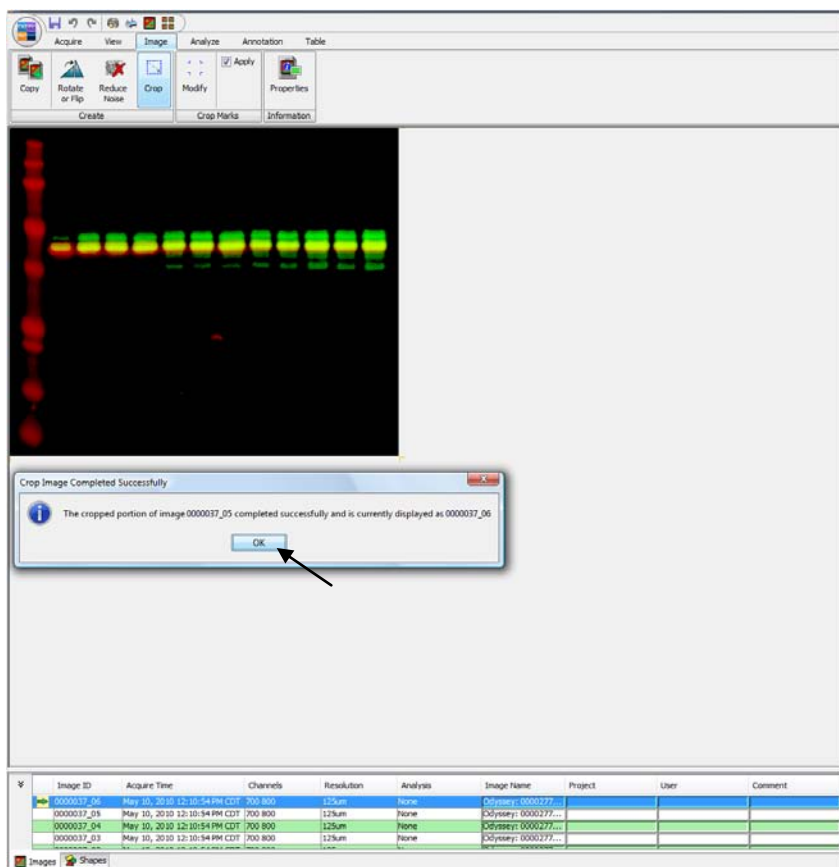


# Chapter 6: Analyzing a Blot or Gel

## Adding and Manipulating Shapes Manually


- 1) Click on the *Analyze* tab to open the *Analyze* ribbon.

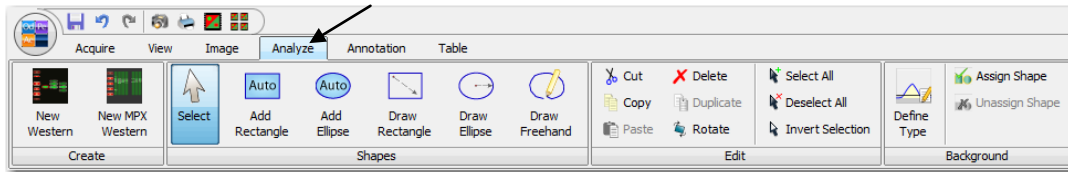


- 2) To manually add shapes to the image, only one channel can be selected. Click on the *Don't show this channel* thumbnail in the Image LUTs for the 700 channel, leaving only the 800 channel on the image.

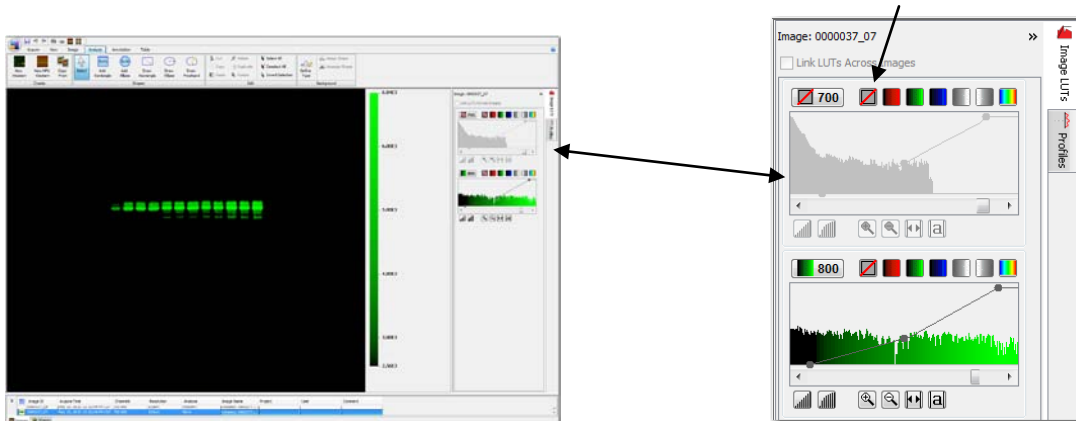

- 3) Click on **Auto Add Rectangle** to add a rectangle to the image or **Auto Add Ellipse** to add an ellipse. Click on the image to place the rectangle or ellipse around a band. After adding the desired shapes, click on **Select** or press the *Esc* key on your keyboard to return the cursor to the selection tool.

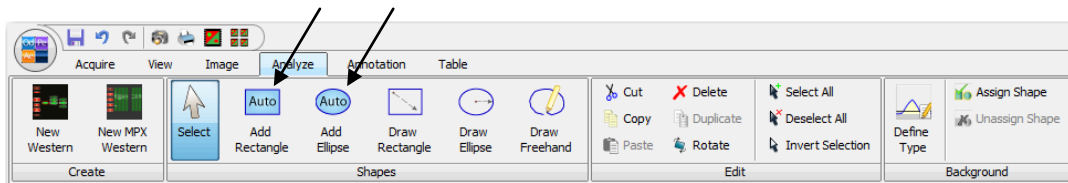


- 4) To move or resize a shape, first select the shape by clicking on it once. Dotted lines indicate the shape is selected.
  - a. To move the shape, click and drag the four-pointed cursor that appears on the shape.

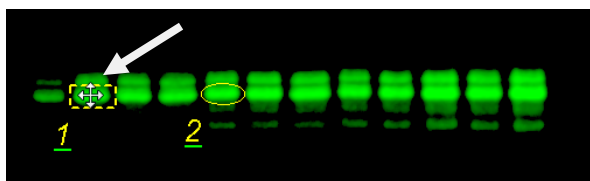

- b. To resize the shape, hover with the mouse over any corner or side of the shape. Click and drag the double-sided yellow arrow to make the shape bigger or smaller.

**Note:** To find the double-sided yellow arrow, the cursor must be on the selection tool (click on **Select** in the Shapes group or push the Esc key on your keyboard).

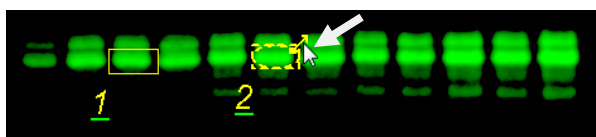

- c. To view the shapes and labels on the image, click on **Shapes** and **Labels** in the View ribbon Show group so that check marks appear in the boxes next to them. Or, remove the check marks to remove them from the image. Refer to [Chapter 5: Manipulating an Image](#).
- d. To edit the label, click on the **Shapes** tab at the bottom of the screen to view the Shapes Table. Under the column **Name**, double-click on the cell for this shape and edit the text. Press **Enter** on the keyboard and the new text will appear on the image. To change the location of the label, refer to [Chapter 5: Manipulating an Image](#).

Image ID	Image Name	Channel	Name	Signal	%
77_01 0000277_01	77_01 0000277_01	800	2	1.13E6	
77_01 0000277_01	77_01 0000277_01	800	2	1.53E6	
77_01 0000277_01	77_01 0000277_01	800	2	1.25E6	
77_01 0000277_01	77_01 0000277_01	800	Text	1.42E6	
77_01 0000277_01	77_01 0000277_01	800	1	1.38E6	
77_01 0000277_01	77_01 0000277_01	700	13	2.00E7	


**Note:** Shapes can be manually drawn by selecting **Draw Rectangle**, **Draw Ellipse**, or **Draw Freehand**. Click and drag the mouse to create the shape. The resulting shape can be moved or resized in the same ways as an auto shape.

- 5) To copy the same shapes to the 700 Channel,

- a. Click **Select All** in the *Edit* group to select the shapes. Dotted lines indicate the shapes are selected.
- b. Click on **Duplicate** in the *Edit* group or right-click on the image and select **Duplicate** from the right-click menu.
- c. Click on the *Show this Channel in Red* thumbnail for the 700 channel to see the bands.



- 6) To copy shapes in the 800 channel...
  - a. Click on the *Don't show this channel* thumbnail for the 700 channel, leaving only the 800 channel on the image.
  - b. Click and drag a rectangle around the shapes to select them. Dotted lines indicate the shapes are selected.
  - c. Click on **Copy** in the *Edit* group or right-click on the image and select **Copy** from the right-click menu.
  - d. Click on the image and click on **Paste** in the *Edit* group or right-click on the image and select **Paste** from the right-click menu.
  - e. To move the shapes, click and drag the four-pointed cursor that appears on either shape. The selected shapes move together.

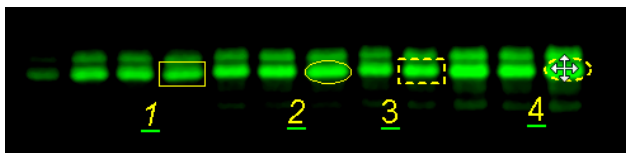


- 7) To rotate the shapes in the 800 channel...
  - a. Click on the *Don't show this channel* thumbnail for the 700 channel, leaving only the 800 channel on the image.
  - b. Click and drag a rectangle around the shapes to select them. Dotted lines indicate the shapes are selected.
  - c. Click on **Rotate** in the *Edit* group or right-click on the image and select **Rotate** from the right-click menu.
  - d. Click and drag the mouse in the circle to rotate the selected shapes.

**Note:** Shapes can also be rotated by selecting them and using the mouse scroll.

## Using the Profiles

- 1) Select 'Shape 2' by clicking on the shape with the selection arrow until it is a dashed line.
- 2) On the right side of the screen, click on the **Profiles** tab to view two graphs of the pixels in a selected shape. The top graph is a cross section of the pixel intensities from top to bottom. The bottom graph is a cross section of the pixel intensities from left to right.

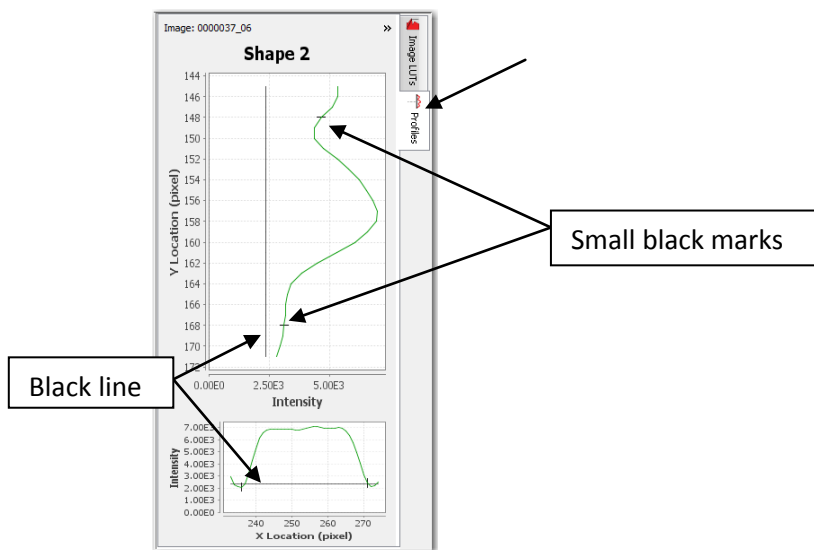


- 3) Using what you've learned from *Adding and Manipulating Shapes Manually* in this chapter, adjust the size or location of 'Shape 2'. Notice how the graph changes.

**Note:** The small black marks indicate the boundaries of the shape. Move the shape around the band so that the small black marks are close to the edges of the peak to capture the entire signal for that band.

## Subtracting the Background

To obtain consistent data, subtract the background of the gel or blot from the shapes.

- 1) Click on **Define Type** in the *Background* group to view the *Background Method* menu.

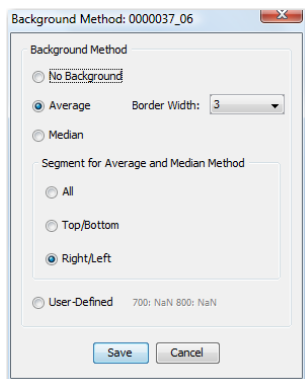



- 2) Click on **Average** to subtract the average value of the pixels in the background segment. Set the *Border Width* to **3**.

**Note:** Select **Median** for the Background Method to subtract the median value of the pixels in the background segment. This method may be useful if the pixel intensities in the background segment are not uniform.


- 3) Click on **Right/Left** to select the segments on the left and right sides of the shape. Other options include selecting the top and bottom sides of the shape or all sides of the shape.
- 4) Click on **Save** to use these settings. A confirmation message will appear if the background method is changed. Click on **Yes** to change the settings.

**Note:** The black line in the **Profiles** indicates the background that will be subtracted.

### Subtracting a User-Defined Background


A region of the image can be chosen for the background.

- 1) Using what you've learned from *Adding and Manipulating Shapes Manually* in this chapter, place a shape on the image in the region that you choose as the background.
- 2) Select the shape by clicking on it with the selection arrow until it is a dashed line.
- 3) Click on **Assign Shape** in the *Background* group. A confirmation message will appear if the background method is changed. Click on **Yes** to change the settings.

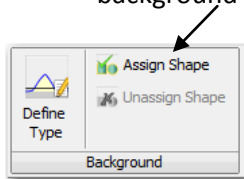


### Annotating the Image


- 1) Click on the **Annotation** tab to open the *Annotation* ribbon.

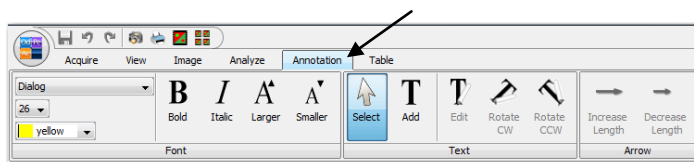

- 2) Click on **Add** to change the cursor to the white text tool. Click anywhere on the image to add text or an arrow at that location. The *Text Properties* menu appears.
- 3) Click on **Text** or **Text with border** in the *Text Properties* menu to add the text inserted in the *Text:* box to the image. Click on **Arrow** to add an arrow to the image.

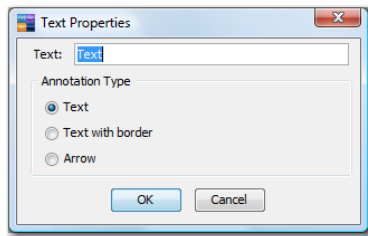

- 4) Click on **Select** or press the *Esc* key on the keyboard to return the cursor to the selection tool.
- 5) Click on the text or arrow that was added to the image with the selection tool. A dashed box appears around the text or the arrow becomes a dashed line.
- 6) Use the tools in the *Font* group to change the look of the text or arrow.

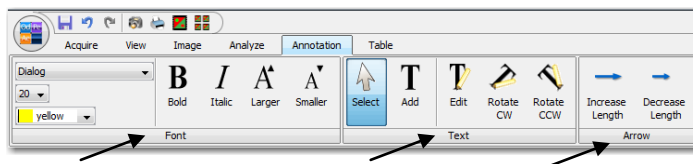

- 7) Click on **Rotate CW** or **Rotate CCW** in the *Text* group to turn the text or arrow.
- 8) Click on **Edit** in the *Text* group to view the *Text Properties* menu for the selected item. Insert a different text or change the text to an arrow in this menu.
- 9) If the selected item is an arrow, click on **Increase** or **Decrease Length** in the *Arrow* group to change the length of the arrow. Change the Font size to change the size of the arrow.


# Chapter 7: Performing Western and MPX™ Western Analyses

---


*Install the Image Studio Western Analysis Key to access the Western Analysis and MPX Western Analysis ribbons. These ribbons have added features for quickly analyzing Westerns and MPX Westerns.*

## Applying a Western Analysis

- 1) Return to the adjusted Western image from [Chapter 5: Manipulating an Image](#) by clicking on the appropriate row in the *Images Table*. To quickly find this image, hover the mouse over the Images Table to view a small picture of the image contained in each row.

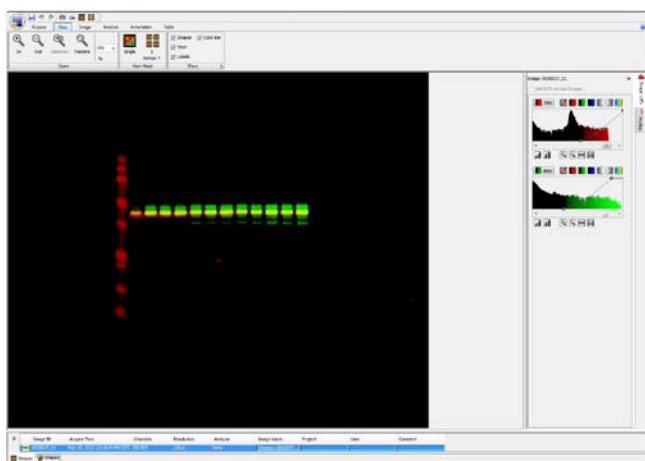


- 2) Click on the **Analyze** tab to view the *Analyze* ribbon.

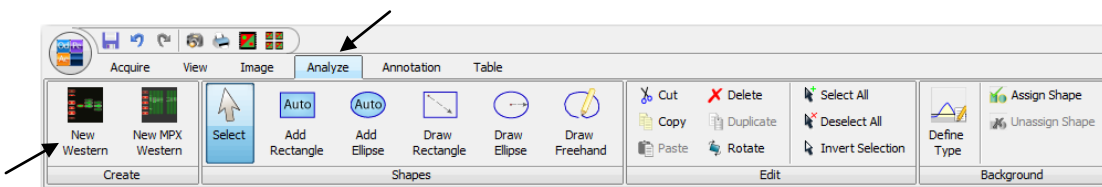

- 3) Click on **New Western** to access the *Western Analysis* ribbon and apply the Western analysis to the image.

**Note:** To apply the Western Analysis automatically after image acquisition, click on **Western** in the Analyze Type group on the Acquire ribbon when starting the acquisition.

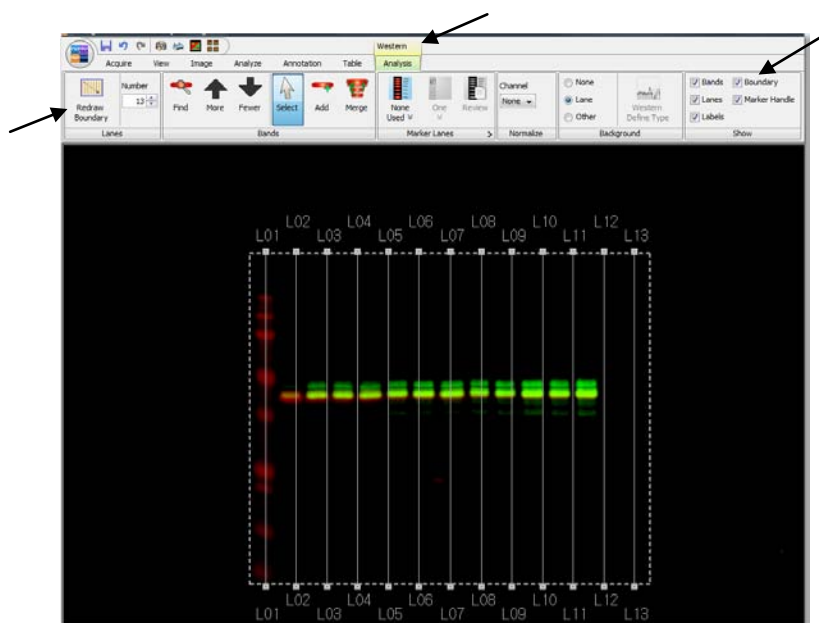

*Settings from the last Western analysis in your Work Area will be applied to the image.*

## Setting the Boundary

*The Boundary is the outside box that surrounds the lanes and bands on the image.*

- 1) To view the Boundary on the image, click on **Boundary** in the **Show** group so that a check mark appears in the box next to it. To hide the Boundary, click to remove the check mark.
- 2) To select the Boundary, click on it once. The line is now dashed indicating it can be adjusted.
- 3) To adjust the applied boundary, hover over the corners or the midpoints of the sides to view a double-sided arrow. Click and drag the arrow to move the corner or side.
- 4) To move the entire Boundary, click and drag the four-pointed cursor that appears anywhere inside the Boundary.
- 5) To draw a new Boundary, click on **Redraw Boundary** in the **Lanes** group. Click and drag a rectangle shape on the image. Repeat the above steps to adjust the Boundary.
- 6) To set the Boundary, click on a different spot on the screen. The Boundary line is now solid.

## Setting the Lanes


*The Lanes are vertical lines that transect the bands in a lane.*

- 1) To view the Lanes on the image, click on **Lanes** in the **Show** group so that a check mark appears in the box next to it. To hide the Lanes, click to remove the check mark.
- 2) Count the total number of lanes in the image (including empty lanes). Adjust the **Number** in the **Lanes** group to match the total number of lanes in the image by clicking on the up and down arrows next to the box.

**Note:** The Lanes are labeled with 'L' plus a number, starting with '01' on the left. To view the labels, click on **Labels** in the Show group so that a check mark appears in the box next to it. To hide the labels, click to remove the check mark.


- 3) To easily adjust the Lanes to transect the image lanes, use the steps in *Setting the Boundary* to adjust the sides of the Boundary.

**Note:** The Lanes are anchored to the Boundary at the top and bottom of the Lane. Adjusting the Boundary also adjusts the Lanes.

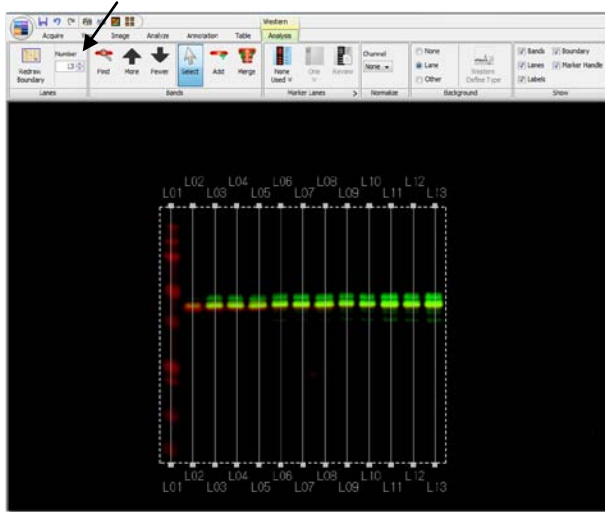

- 4) To further adjust a single Lane, click on it once so that the line appears dashed. Click and drag the small solid boxes at the top and bottom of the Lane along the Boundary.

**Note:** The Lanes only need to be close to the centers of the bands, as they will be optimally adjusted by the software during band finding.

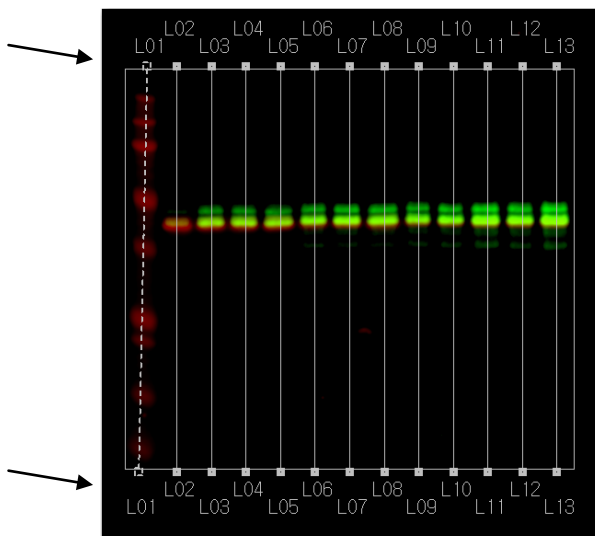

- 5) To set the Lane, click on a different spot on the screen. The Lane line is now solid.


## Setting the Marker

*Each Marker Lane is designated by 'M' followed by 1, 2, or 3.*

- 1) Click on the *Select Marker Lanes Set* icon in the *Marker Lanes* group and choose **One-Color Protein** from the pull-down menu.
- 2) Click on the *Number of Marker Lanes* icon in the *Marker Lanes* group and select **One** from the pull-down menu.
- 3) To change the location of the Marker Lane, click on the Marker Handle to change the rectangle to a dashed line.

**Note:** *The Marker Handle is a rectangle with 'M1', 'M2', or 'M3' inside. It designates the lane adjacent to its left side as a marker lane. To view the Marker Handle, click on **Marker Handle** in the Show group so that a check mark appears in the box next to it. To hide the Marker Handle, click to remove the check mark.*

- 4) Click and drag the four-pointed cursor on the Marker Handle over any lane to make that lane the Marker. For the tutorial, drag the cursor to the first lane to make it the Marker. The name of the lane changes to 'M1' and the other lanes are renumbered to exclude that lane.

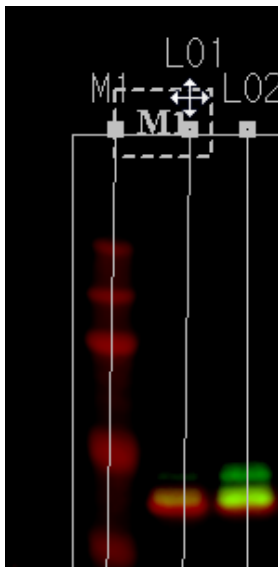

- 5) Click on **Review** in the *Marker Lanes* group to see the assigned values for the bands in the One-Color Protein Marker. Click **OK** to close the box.

## Creating a New Marker

- 1) Click on the expand button on the *Marker Lanes* group to open the *Create New or Delete Marker* menu.

- 2) Click on **New...** to open the *Name Marker* menu. Enter a name for the Marker in the box and click **OK**. The *New Marker* menu will open.
- 3) Click on **Add** to open the *Add Band Value* menu. Enter a numerical value for a Marker band in the box and click **OK**.
- 4) Repeat step 3 to add all of the values for the bands in the Marker. The numerical values will be sorted automatically from high to low. To delete a value, click on the value to select it and click on **Delete**.
- 5) Click on the down arrow next to *Units* to view the pull-down menu. Click on the desired units from the options shown.
- 6) In the Channels group, click on **Add** to view the Channel Name menu. Enter 700, 800, or Chemi and click **OK**. Repeat to enter multiple channels. To delete an entry, click on the entry to select it and click on **Delete**.
- 7) Click **OK** to close the *New Marker* menu.

**Note:** The name of the new Marker will appear in the Select Marker Lanes Set pull-down menu in the Marker Lanes group. Select the Marker from the menu and click on **Review** to see the details. Click on the expand button on the Marker Lanes group to open the Create New or Delete Marker menu and click on **Delete** to remove the Marker.


## Finding the Bands Automatically

*Bands will be found in all channels.*

- 1) In the *Bands* group, click on **Find** to automatically find bands on the Lanes.

**Note:** Removing the Labels makes the Bands easier to see. Click on **Labels** in the Show group to remove the check mark next to it.

- 2) Click on **Fewer** to decrease the sensitivity resulting in fewer Bands. Click on **More** to increase the sensitivity and find more Bands.

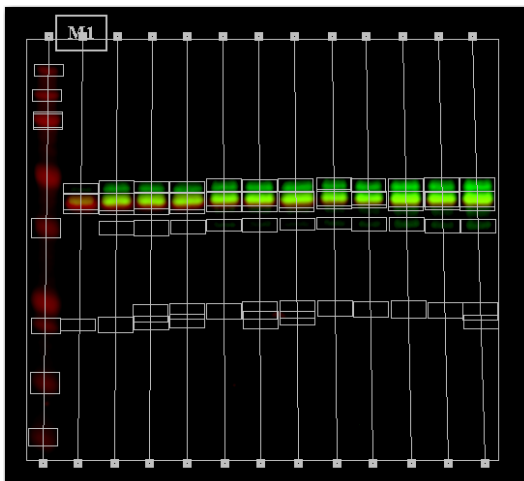


## Editing Bands Manually

*Only one channel may be selected to edit Bands manually.*

- 1) To manually edit the Bands, only one channel can be selected. Click on the **Don't show this channel** thumbnail for the 800 channel, leaving only the 700 channel on the image.
- 2) To manually add a Band,
  - a. Click once on **Add** in the *Bands* group and then click once on a Lane to place the Band.
  - b. Click on **Select** or press the *Esc* key on the keyboard to return the cursor to the selection tool.
- 3) To manually delete a Band,
  - a. Select the Band by clicking on it once with the selection tool. Dotted lines indicate the Band is selected.
  - b. Press *Ctrl+X* or the *Delete* key on the keyboard (or right-click and select **Cut** or **Delete**) to delete the selected Band.

**Note:** Multiple Bands can be deleted or moved at one time. Select multiple Bands by pressing the *Ctrl* key on the keyboard and clicking once on each Band or use the selection tool to click and drag a rectangle over the Bands.

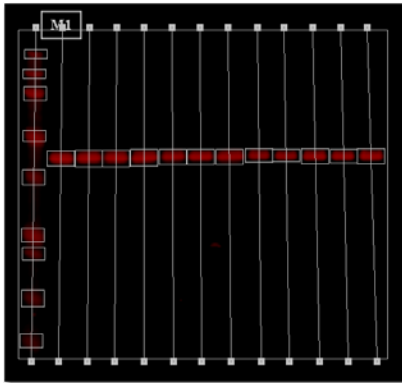

- 4) To move a Band,
  - a. Select the Band by clicking on it once with the selection tool. Dotted lines indicate the Band is selected.
  - b. Click and drag the four-pointed cursor that appears on the selected Band.

**Note:** The Band is anchored to the Lane it appears on and cannot be moved to another Lane.

- 5) To resize a Band,
  - a. Select the Band by clicking on it once with the selection tool. Dotted lines indicate the Band is selected.
  - b. Hover with the mouse over any side of the Band. Click and drag the double-sided arrow to make the selected Band bigger or smaller.
- 6) To merge Bands,



- a. Use the selection tool to click and drag a rectangle over two or more Bands that appear on the same Lane, or press *Ctrl* on the keyboard and click once on each Band. Dotted lines indicate the Bands are selected.
- b. Click on **Merge** in the *Band* group. One Band now surrounds the area in and between the previous Bands.

## Subtracting the Background

To obtain consistent data, subtract the background of the blot from the Bands.

- 1) Click on **Lane** in the *Background* group to subtract the mean background in the Lane from the Bands that appear on that Lane.

**Note:** *Lane* is the recommended setting for background subtraction in the Western Analysis application.

- 2) To define a different background subtraction, click on **Other** in the *Background* group to enable the **Western Define Type** icon.
- 3) Click on **Western Define Type** to view the *Western Background* menu. Refer to *Subtracting the Background* in [Chapter 6: Analyzing a Blot or Gel](#) for more information on this menu.

## Reviewing the Lanes Table

- 1) Click on the **Lanes** tab at the bottom of the screen to view the Lanes Table. The Lanes Table shows the Lanes from the current single image.

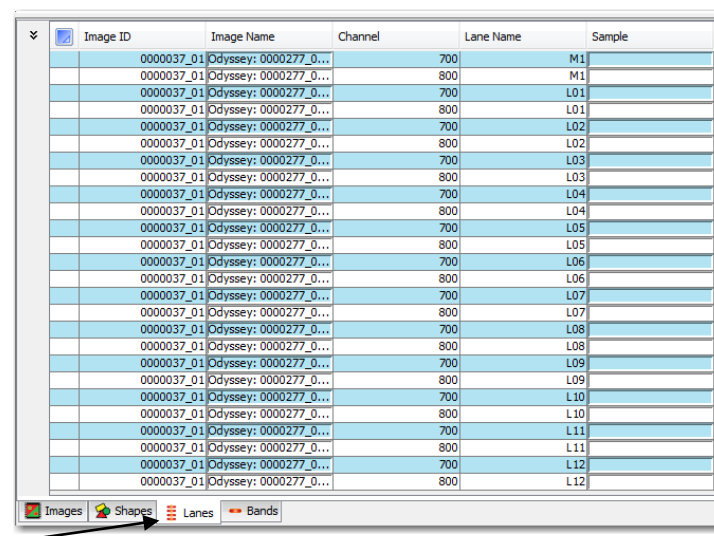


Image ID	Image Name	Channel	Lane Name	Sample
0000037_01	Odyssey: 0000277_0...	700	M1	
0000037_01	Odyssey: 0000277_0...	800	M1	
0000037_01	Odyssey: 0000277_0...	700	L01	
0000037_01	Odyssey: 0000277_0...	800	L01	
0000037_01	Odyssey: 0000277_0...	700	L02	
0000037_01	Odyssey: 0000277_0...	800	L02	
0000037_01	Odyssey: 0000277_0...	700	L03	
0000037_01	Odyssey: 0000277_0...	800	L03	
0000037_01	Odyssey: 0000277_0...	700	L04	
0000037_01	Odyssey: 0000277_0...	800	L04	
0000037_01	Odyssey: 0000277_0...	700	L05	
0000037_01	Odyssey: 0000277_0...	800	L05	
0000037_01	Odyssey: 0000277_0...	700	L06	
0000037_01	Odyssey: 0000277_0...	800	L06	
0000037_01	Odyssey: 0000277_0...	700	L07	
0000037_01	Odyssey: 0000277_0...	800	L07	
0000037_01	Odyssey: 0000277_0...	700	L08	
0000037_01	Odyssey: 0000277_0...	800	L08	
0000037_01	Odyssey: 0000277_0...	700	L09	
0000037_01	Odyssey: 0000277_0...	800	L09	
0000037_01	Odyssey: 0000277_0...	700	L10	
0000037_01	Odyssey: 0000277_0...	800	L10	
0000037_01	Odyssey: 0000277_0...	700	L11	
0000037_01	Odyssey: 0000277_0...	800	L11	
0000037_01	Odyssey: 0000277_0...	700	L12	
0000037_01	Odyssey: 0000277_0...	800	L12	

**Note:** The columns in the Lanes Table can be moved, added, deleted, or sorted in the same ways as the columns in the Images Table. The Lanes Table cannot be filtered. Refer to [Chapter 8: Working with the Tables](#) for more information.

## Reviewing the Bands Table

- 1) Click on the **Bands** tab at the bottom of the screen to view the Bands Table. The Bands Table shows the Bands from the current single image.


Image ID	Image Name	Channel	Lane Name	Name	Signal	Size
0000037_01	Odyssey: 0000277_0...	700	M1	B01	6.24E5	2.50E2
0000037_01	Odyssey: 0000277_0...	700	M1	B02	8.50E5	1.50E2
0000037_01	Odyssey: 0000277_0...	700	M1	B03	1.44E6	1.00E2
0000037_01	Odyssey: 0000277_0...	700	M1	B04	1.18E6	7.50E1
0000037_01	Odyssey: 0000277_0...	700	M1	B05	8.81E5	5.00E1
0000037_01	Odyssey: 0000277_0...	700	M1	B06	1.69E6	3.70E1
0000037_01	Odyssey: 0000277_0...	700	M1	B07	7.60E5	2.59E1
0000037_01	Odyssey: 0000277_0...	700	M1	B08	1.02E6	2.00E1
0000037_01	Odyssey: 0000277_0...	700	M1	B09	9.10E5	1.50E1
0000037_01	Odyssey: 0000277_0...	800	M1	B01	1.18E5	2.50E2
0000037_01	Odyssey: 0000277_0...	700	L01	B01	3.00E6	6.02E1
0000037_01	Odyssey: 0000277_0...	800	L01	B01	2.61E5	6.83E1
0000037_01	Odyssey: 0000277_0...	800	L01	B02	6.74E5	6.12E1
0000037_01	Odyssey: 0000277_0...	700	L02	B01	3.20E6	6.02E1
0000037_01	Odyssey: 0000277_0...	800	L02	B01	7.74E5	6.89E1
0000037_01	Odyssey: 0000277_0...	800	L02	B02	1.49E6	6.12E1
0000037_01	Odyssey: 0000277_0...	800	L02	B03	7.10E4	5.00E1
0000037_01	Odyssey: 0000277_0...	700	L03	B01	3.21E6	6.02E1
0000037_01	Odyssey: 0000277_0...	800	L03	B01	7.13E5	6.89E1
0000037_01	Odyssey: 0000277_0...	800	L03	B02	1.54E6	6.17E1
0000037_01	Odyssey: 0000277_0...	800	L03	B03	9.04E4	5.00E1
0000037_01	Odyssey: 0000277_0...	800	L03	B04	1.09E5	3.13E1
0000037_01	Odyssey: 0000277_0...	700	L04	B01	3.42E6	6.07E1
0000037_01	Odyssey: 0000277_0...	800	L04	B01	8.22E5	6.89E1
0000037_01	Odyssey: 0000277_0...	800	L04	B02	1.59E6	6.12E1
0000037_01	Odyssey: 0000277_0...	800	L04	B03	7.89E4	5.03E1

**Note:** The columns in the Bands Table can be moved, added, deleted, or sorted in the same ways as the columns in the Images Table. The Bands Table cannot be filtered. Refer to [Chapter 8: Working with the Tables](#) for more information.

- 2) Notice the first 10 Bands are in the Marker Lane (M1) and the values from the selected Marker are in the column Size. The rest of the Bands are assigned a Size in reference to the values of the Marker Bands.


## Applying the MPX™ Western Analysis

- 1) Import 'MPX data.zip' from the Image Examples folder on the Image Studio software CD. Refer to [Chapter 4: Importing an Image](#) for more information.
  - a. Click on the **Application Menu** button and move the mouse to hover over **Import**. Click on **Copy From (Import Acquisition)** in the **Import** menu.
  - b. Click once on the file '9999998\_01' to select it. Click **Open** to display the image. The file will appear in the Images Table.
- 2) Adjust the look of the table in the **Image LUTs** on the right side of the screen. Refer to [Chapter 5: Manipulating an Image](#) for more information.
- 3) Click on the **Analyze** tab to view the **Analyze** ribbon.

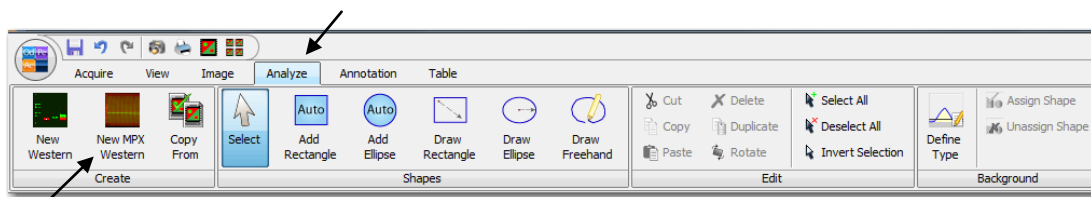


- 4) Click on **New MPX Western** in the **Create** group to open the **MPX Western Analysis** ribbon and apply the MPX Western analysis to the image.

**Note:** To apply the MPX™ Western analysis automatically after image acquisition, click on **MPX Western** in the Analyze Type group on the Acquire ribbon when starting the acquisition.

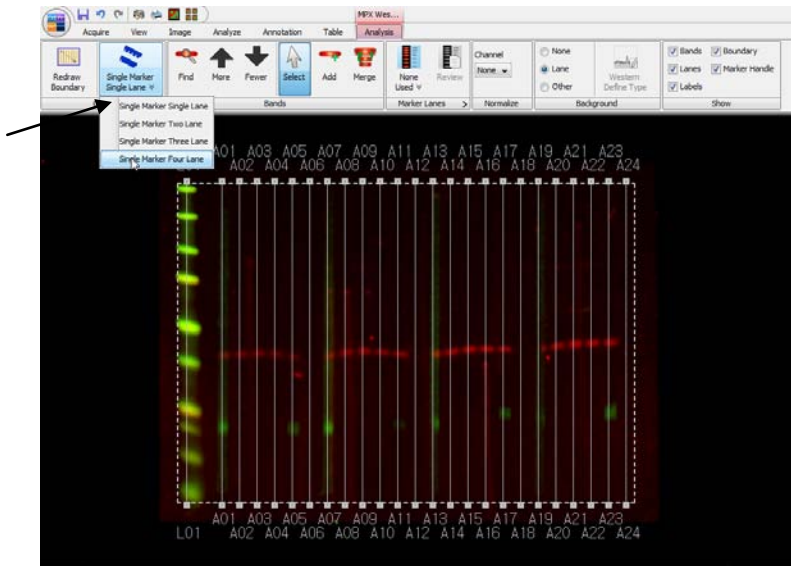

Parameters from the last MPX Western analysis will be applied to the image.

- 5) Click on **Select Comb for MPX Western** to open the pull-down menu. Click on **Single Marker Four Lane** to apply a template with a single marker followed by four sets of five Lanes each.

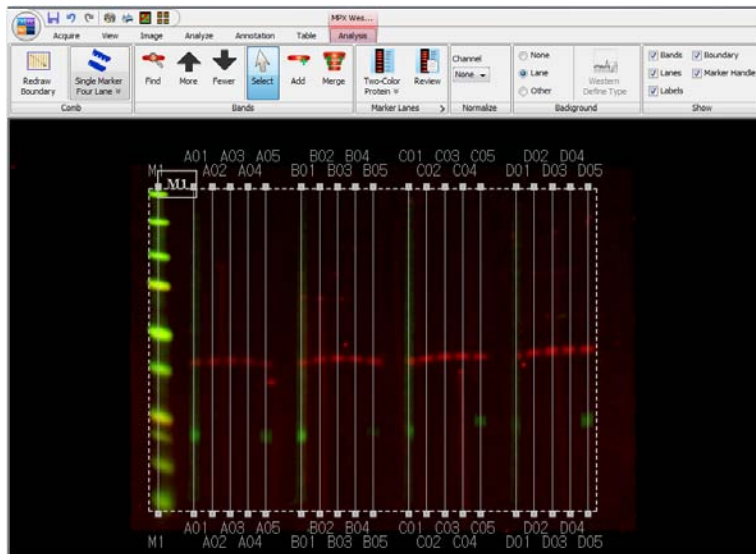

- 6) Adjust the Boundary so the Lanes transect the image lanes. Refer to *Setting the Boundary* in this chapter for more information.
- 7) Click on **Select Marker Lanes Set** in the *Marker Lanes* group to view the pull-down menu. Click on **Two-Color Protein**. Refer to *Setting the Marker* in this chapter for more information.

- 8) Click on **Find** in the *Bands* group to find the Bands automatically. Refer to *Finding the Bands Automatically* and *Editing Bands Manually* in this chapter for more information.
- 9) Click on **Lane** in the *Background* group to set the background subtraction. Refer to *Subtracting the Background* in this chapter for more information.

## Reviewing the MPX™ Lanes Table

- 1) Click on MPX Lanes at the bottom of the screen to view the MPX Lanes Table.

Image ID	Image Name	Channel	Lane Name	Sample
0000418_01	0000418_01	7960	M1	
0000418_01	0000418_01	8000	M2	
0000418_01	0000418_01	7960	M3	
0000418_01	0000418_01	8000	M4	
0000418_01	0000418_01	7960	M5	
0000418_01	0000418_01	8000	M6	
0000418_01	0000418_01	7960	M7	
0000418_01	0000418_01	8000	M8	
0000418_01	0000418_01	7960	M9	
0000418_01	0000418_01	8000	M10	
0000418_01	0000418_01	7960	M11	
0000418_01	0000418_01	8000	M12	
0000418_01	0000418_01	7960	M13	
0000418_01	0000418_01	8000	M14	
0000418_01	0000418_01	7960	M15	
0000418_01	0000418_01	8000	M16	
0000418_01	0000418_01	7960	M17	
0000418_01	0000418_01	8000	M18	
0000418_01	0000418_01	7960	M19	
0000418_01	0000418_01	8000	M20	
0000418_01	0000418_01	7960	M21	
0000418_01	0000418_01	8000	M22	


**Note:** The columns in the MPX Lanes Table can be moved, added, deleted, or sorted in the same ways as the columns in the Images Table. The MPX Lanes Table cannot be filtered. Refer to [Chapter 8: Working with the Tables](#) for more information.

# Chapter 8: Working with the Tables


Image data is stored in the folder selected as the Work Area when the Image Studio software is opened. The data is easily accessed in the Tables below the image window. In the Images Table, each row represents one image. Various fields in that row contain identifying information about the image. These fields are grouped as columns that can be moved, added, deleted, sorted, or filtered.

## Working in the Images Table

- 1) Click on the **Images** tab at the bottom of the screen to show the Images Table. (Click on the Expand button to hide the Table from view.)

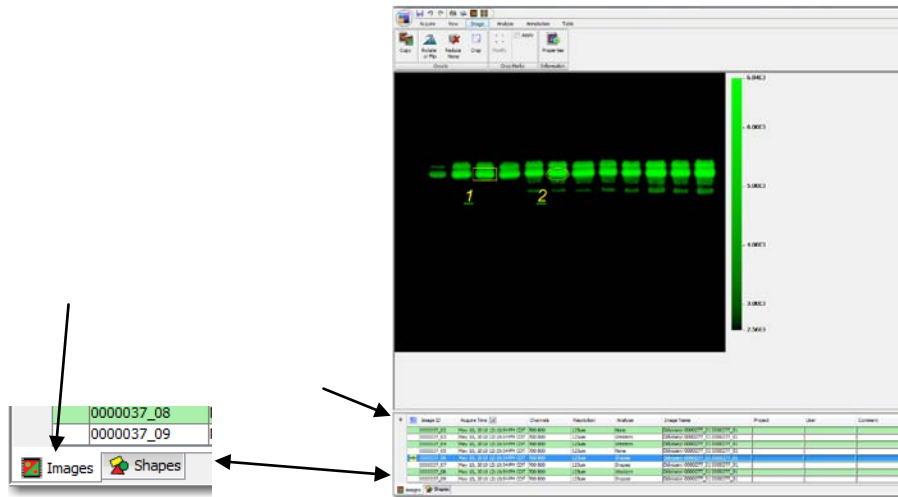


- 2) Hover with the mouse over any row to show a thumbnail picture of that image.

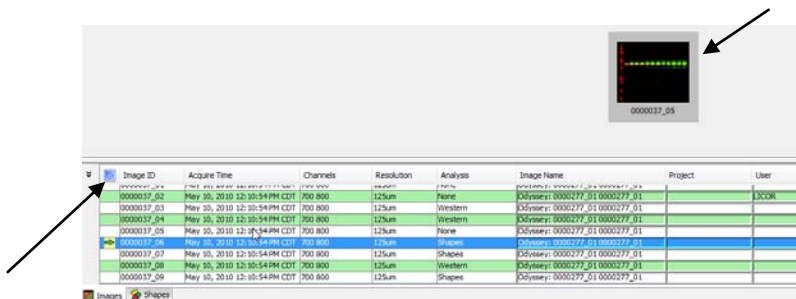

- 3) Click on any cell to select the entire row and show that image in the window.
- 4) Click on the corner button to select all of the images or press **Ctrl+A** on the keyboard.
- 5) Right-click any cell to view a menu with options that include copy an image, delete images, or view the properties of an image.

- 6) To move a column, click and drag the column header to the new location.

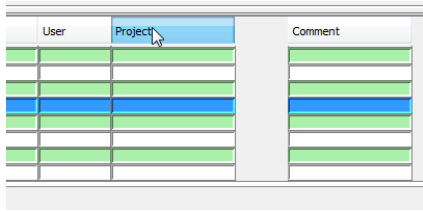


*The fields in each row contain information about the image and are organized as columns across all the images. Some of the columns are properties of the image and cannot be edited. Examples of these columns include Image ID, Acquire Time, Channels, Resolution, and Instrument Name. Other columns can be edited with information for that image.*

- 7) To edit a field, double-click on a cell so that a cursor appears. If a cursor does not appear, that field cannot be edited.
- 8) Add text where the cursor appears. Press *Enter* on the keyboard when complete.


**Note:** To easily find images later, enter identifying information in one or more fields after acquiring an image.

## Adding or Deleting a Column

- 1) Click on the **Table** tab to view the *Table* ribbon.
- 2) Click on **Add/Remove** in the *Columns* group.

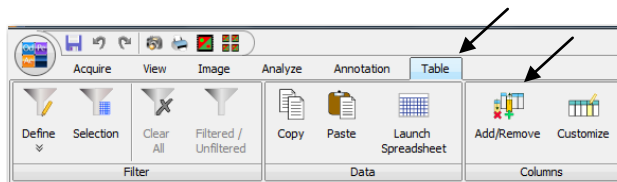

- 3) To add the *Experiment* column, click the column description to show a check mark indicating that column will be in the table.
- 4) To remove the *Resolution* column, click the column description to remove the check mark and remove that column from the table.

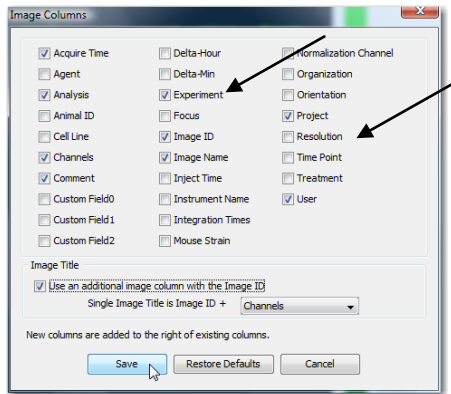


5) Click on **Save** to use the changes. The table now looks like this one.


Image ID	Acquire Time	Channels	Analysis	Image Name	User	Project	Comment	Experiment
0000037_06	May 10, 2010 12:10:54 PM CDT	700 800	Shapes	Odyssey: 0000277_01 0000277_01				
0000037_07	May 10, 2010 12:10:54 PM CDT	700 800	Shapes	Odyssey: 0000277_01 0000277_01				
0000037_09	May 10, 2010 12:10:54 PM CDT	700 800	Shapes	Odyssey: 0000277_01 0000277_01				
0000037_01	May 10, 2010 12:10:54 PM CDT	700 800	None	Odyssey: 0000277_01 0000277_01				
0000037_02	May 10, 2010 12:10:54 PM CDT	700 800	None	Odyssey: 0000277_01 0000277_01				
0000037_05	May 10, 2010 12:10:54 PM CDT	700 800	None	Odyssey: 0000277_01 0000277_01				
0000038_01	May 14, 2010 9:11:07 AM CDT	700 800	None	Odyssey: 0000002_01 0000002_01				
0000039_01	May 14, 2010 3:10:33 PM CDT	700 800	None	Odyssey: 0000002_01 0000002_01				

## Editing the Image Title

- 1) Click on the **Table** tab to view the *Table* ribbon.
- 2) Click on **Add/Remove** in the *Columns* group.

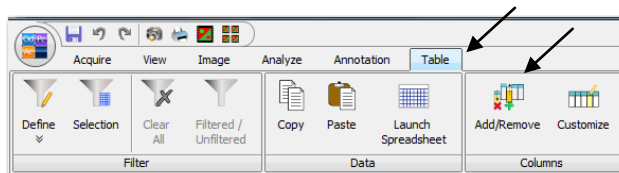


- 3) Click on *Use an additional image column with the Image ID* to place a check mark in the check box next to it.

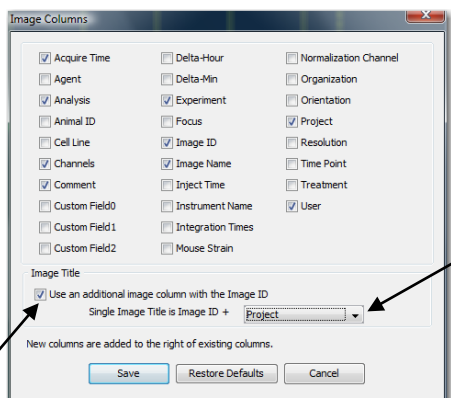


- 4) Click on the down arrow in the selection box to choose a column to add to the Image ID to form the Image Title.
- 5) Click on **Project**. Text that is in the *Project* column will be added to the title of every Image.
- 6) Click on **Save** to use the changes.

## Adding Custom Columns


- 1) Click on the **Table** tab to view the *Table* ribbon.
- 2) Click on the **Customize** icon in the *Columns* group to view the *Custom Columns* menu.

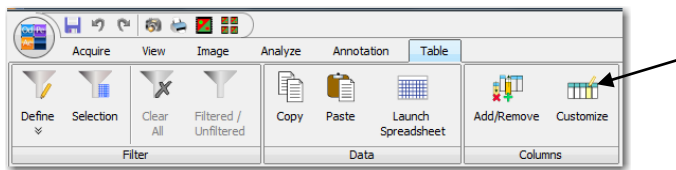

- 3) In the *Custom Columns* menu, replace *Custom Field0* with *Department*.

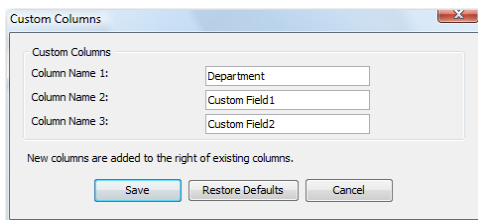

- 4) Click on **Save** to use this change.
- 5) To add this custom column to the table, first click on the **Table** tab to view the *Table* ribbon.
- 6) Click on **Add/Remove** in the *Columns* group.

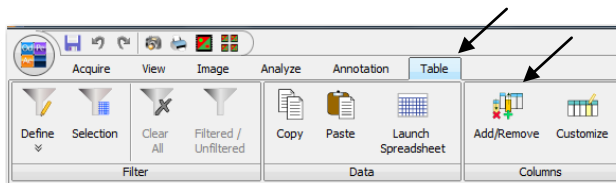

- 7) To add the *Department* column, click the column description to show a check mark indicating that column will be in the table.
- 8) Click on **Save** to use this change. The *Department* column will be added to the right of the existing columns.


## Pasting Data in a Column

Use **Paste** in the *Data* group to quickly fill a column with information.

- 1) Double-click the first cell in the *Department* column so that a cursor appears in the cell.
- 2) Type 'Tutorial' at the cursor position using the keyboard.



- 3) Click and drag the mouse over 'Tutorial' to highlight it. The cursor should still appear in the cell.
- 4) Press **Ctrl+C** on the keyboard to copy the data to the clipboard.
- 5) Click and drag the mouse over the cells in the *Department* column (or press the **Ctrl** key while clicking each cell once) to select which cells to fill with 'Tutorial'.
- 6) Click on **Paste** in the *Data* group, press **Ctrl+V** on the keyboard, or use the right-click menu to paste the data in the selected cells.

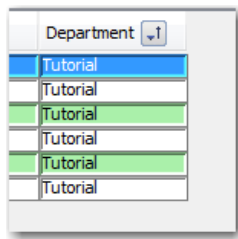

*'Tutorial' added to cells in Department column.*

### Pasting Data from a Spreadsheet to a Column

- 1) In a spreadsheet, select data and press **Ctrl+C** on the keyboard to copy the data to the clipboard.
- 2) In the Table, click and drag the mouse over the cells in the *Department* column (or press the **Ctrl** key while clicking each cell once) to select which cells in a column to place the data.
- 3) Click on **Paste** in the *Data* group, press **Ctrl+V** on the keyboard, or use the right-click menu to paste the data in the selected cells.

### Copying Data to a Spreadsheet

Use **Copy** in the *Data* group to quickly copy rows to a spreadsheet.

- 1) Click and drag the mouse over the rows or press the **Ctrl** key while clicking on each row once to select which rows to copy. To copy all rows, press **Ctrl+A** on the keyboard.
- 2) Click on **Copy** in the *Data* group, press **Ctrl+C** on the keyboard, or use the right-click menu to copy the data to the clipboard.
- 3) In a spreadsheet, press **Ctrl+V** on the keyboard to paste the data in the selected cells.

**Note:** The column headings will be pasted above the data.

### Launching a Spreadsheet

Use **Launch Spreadsheet** in the *Data* group to open an '.xls' or '.xlsx' file with the selected data.

- 1) Click and drag the mouse over the rows or press the **Ctrl** key while clicking on each row once to select which rows to copy.
- 2) Click on **Launch Spreadsheet** to open the *Launch Spreadsheet* browse window.
- 3) Select the folder destination and name the file with extension '.xls' or '.xlsx'. Click **Save**.
- 4) An application that uses these types of files will open with the data selected and corresponding column headings.

## Sorting by Column

- 1) To sort by any column, click once on the header. The table will be sorted by that column.

**Note:** No change will be made if the fields in a column are all the same.

- 2) To change the direction of sorting, click on the small box in the header with the up or down arrow. Or, right-click on the header and select **Sort Ascending** or **Sort Descending**.


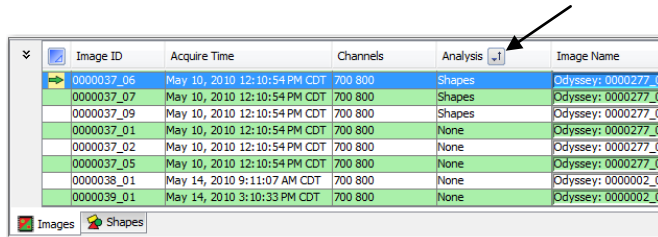



Image ID	Acquire Time	Channels	Analysis	Image Name
0000037_06	May 10, 2010 12:10:54 PM CDT	700 800	Shapes	Odyssey: 0000277_01
0000037_07	May 10, 2010 12:10:54 PM CDT	700 800	Shapes	Odyssey: 0000277_01
0000037_09	May 10, 2010 12:10:54 PM CDT	700 800	Shapes	Odyssey: 0000277_01
0000037_01	May 10, 2010 12:10:54 PM CDT	700 800	None	Odyssey: 0000277_01
0000037_02	May 10, 2010 12:10:54 PM CDT	700 800	None	Odyssey: 0000277_01
0000037_05	May 10, 2010 12:10:54 PM CDT	700 800	None	Odyssey: 0000277_01
0000038_01	May 14, 2010 9:11:07 AM CDT	700 800	None	Odyssey: 0000002_01
0000039_01	May 14, 2010 3:10:33 PM CDT	700 800	None	Odyssey: 0000002_01

## Filtering the Table Manually

- 1) To manually choose images, click and drag on the rows to select the desired images. Or, select the desired images by pressing *Ctrl* on the keyboard and clicking once on each row.
- 2) Click on the **Table** tab to choose the *Table* ribbon.

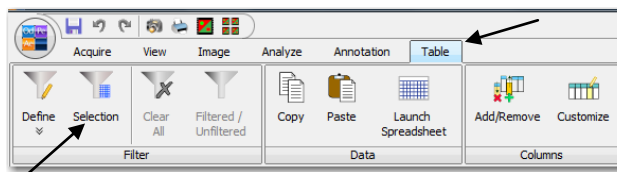

- 3) Click on **Selection** to show only the selected images in the table. The **Filtered/Unfiltered** icon will be highlighted.
- 4) Click on the **Filtered/Unfiltered** icon to toggle between showing only the selected images and showing all images in the table.

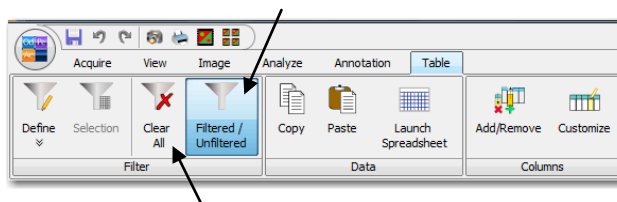

- 5) To remove the filter, click on **Clear All**.


## Defining Filters

One or more defined filters can be applied to the table to show only the images of interest.

- 1) Click on the **Table** tab to choose the *Table* ribbon.
- 2) Click on **Define** to view the pull-down menu of all of the column headings. Select **Analysis** to filter in the *Analysis* column.

**Note:** Right-click on any column header in the Table to access this menu for that column.


- 3) Deselect **None** and **Western** by clicking on the box next to them to remove the check mark.

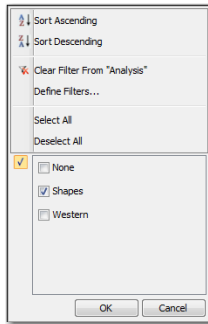

- 4) Click **OK** to use this change. Only the images with *Shapes* in the *Analysis* column will appear in the Table.

### For more Filtering Options...


- 1) Click on **Define Filters...** in the *Filter* menu.
- 2) Click on one of the filtering phrases to show the *Column Filter Analysis* menu.
- 3) Select filtering phrases from the pull-down menu by clicking on the down arrows next to each box or type in your own.

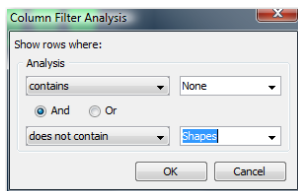

- 4) Click **OK** to use these changes.

## Viewing the Shapes Table

- 1) Click on the **Shapes** tab in the Table to view the data for each shape.

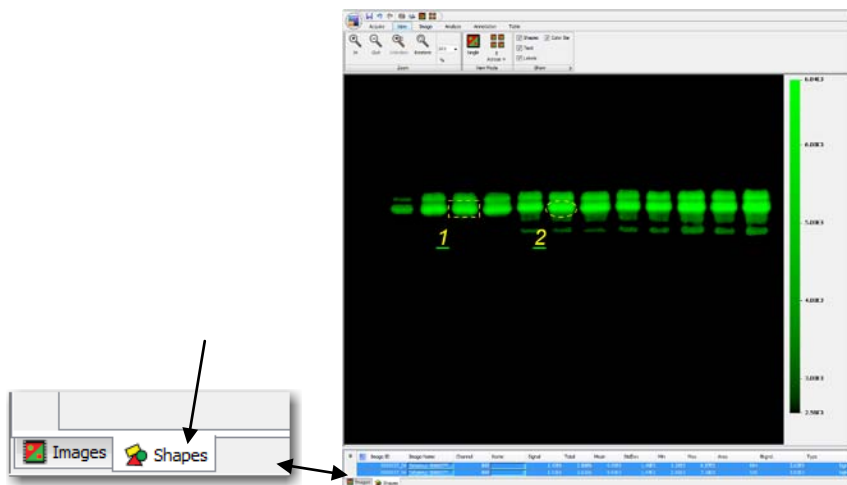

- 2) The default columns are defined below.
  - a. *Image ID*- Number assigned by the software to the image; a new one is assigned whenever a new image is created by cropping, copying, rotating, filtering, or analyzing.
  - b. *Image Name*- Original acquisition name
  - c. *Channel*- 700, 800, or Chemi
  - d. *Name*- Shapes are assigned a number in sequential order. Double-click on the number in the column to edit the shape name.
  - e. *Signal*- Sum of the individual pixel intensities for the shape minus the background.
  - f. *Total*- Sum of the individual pixel intensities for the shape
  - g. *Mean*- Mean of all pixel intensities for a shape
  - h. *StdDev*- Standard deviation of all of the pixel intensities for a shape
  - i. *Min*- Minimum pixel intensity for a shape
  - j. *Max*- Maximum pixel intensity for a shape
  - k. *Area*- Total number of pixels enclosed by a shape
  - l. *Bkgnd*.- Designated background subtraction
  - m. *Type*- Signal or Background
- 3) Move, add, delete, and sort the columns in the same way as in the **Images** Tab. The **Shapes** Tab cannot be filtered.

# Chapter 9: Exporting Data, Printing Images, and Generating Reports

---

## Exporting Acquisition Data and Images

- 1) In the Images Table, click on a row to select that image acquisition. To select multiple image acquisitions, click and drag over the rows or press the *Ctrl* key on the keyboard and click once on each row to select it.
- 2) Click on the **Image Studio Application Button** and hover over **Export** to view the *Export image and analysis data* menu.



- a. To make a copy in another folder, click on **Copy To** and view a choice of **Backup folder** or **Zip File**. Click on **Zip File** to save the image acquisition as a zip file. Either choice opens to a browse window where you can select the folder to copy to.
- b. To make a copy in another folder and delete the image acquisition from the Work Area, click on **Move To (Archive)** and click on **Yes** in the *Move Selected Acquisitions* menu. Select the folder to copy to in the browse window.
- c. To make an image file, click on **Image** to view a choice of **Current Image** or **Selected Images in Images Table**. Either choice opens to a browse window where you can select the folder to copy to as well as the file type (TIFF, PNG, or High Quality JPEG) and the resolution.

**Note:** Images saved as these file types (TIFF, PNG, or High Quality JPEG) are for presentation purposes and should not be used for further analysis.

- d. To make a tiled image file of the images in the Table, click on **Multiple Image View** to open a browse window where you can select the folder to copy to as well as the file type (TIFF, PNG, or High Quality JPEG) and the resolution. In addition, select


**Export Images Table** to copy a Tab Delimited File of the Images Table to the same folder.

- e. To copy a Tab Delimited File of the current Table, click on **Table** to view a choice of **All Rows in Table** or **Selected Rows in Table**. Either choice opens to a browse window where you can select the folder to copy to.


**Note:** Data can be quickly and easily exported from the Tables to an Excel spreadsheet. Refer to Copying Data to a Spreadsheet in Chapter 8: Working with the Tables.

## Printing Images

- 1) To open the *Print* menu, click on the printer icon in the quick access tool bar, press Ctrl+P on the keyboard, or click on the **Image Studio Application Button**, hover over **Print**, and click on **Print**.

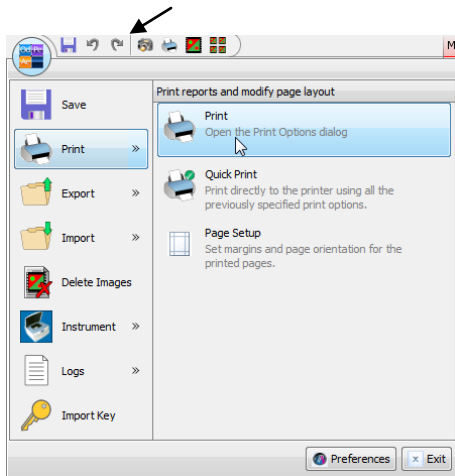


- 2) Click on **Print** to send the current image to your default printer.

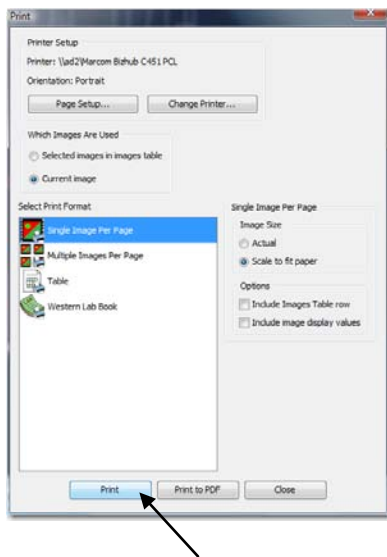


## Generating Reports

- 1) In the *Print* menu, click on **Western Lab Book** and then **Print** to print the default report for the current image.

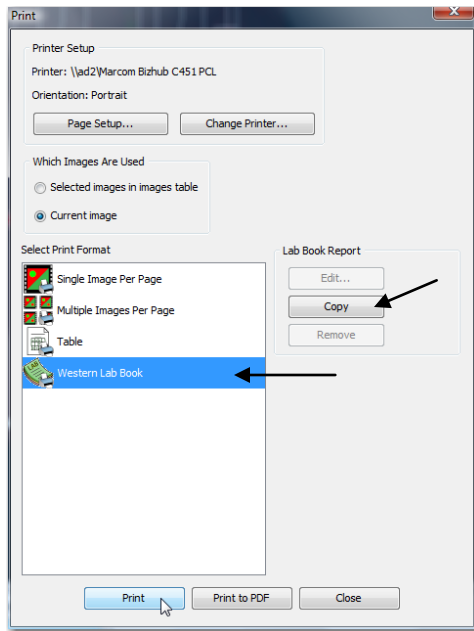


- 2) To create a custom report,
  - a. Click on **Western Lab Book** in the *Print* menu. In the right column under *Lab Book Report*, click on **Copy**.
  - b. In the *Specify Custom Format Print Title* menu, insert 'Tutorial' using the keyboard and click **OK**. **Tutorial** is now a selection in the *Select Print Format* view.

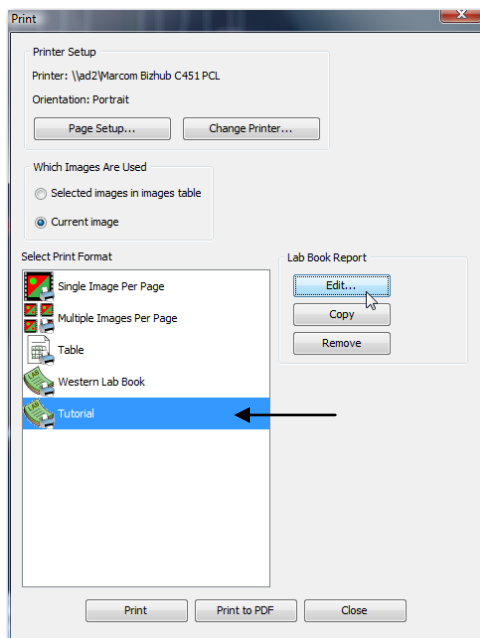

- c. Click on **Edit...** under **Lab Book Report** to see the **Print Format builder** menu.

**Note:** The items currently in the report appear under *In Use* in the left column. Items that are available to add to the report appear under *Available* in the right column.

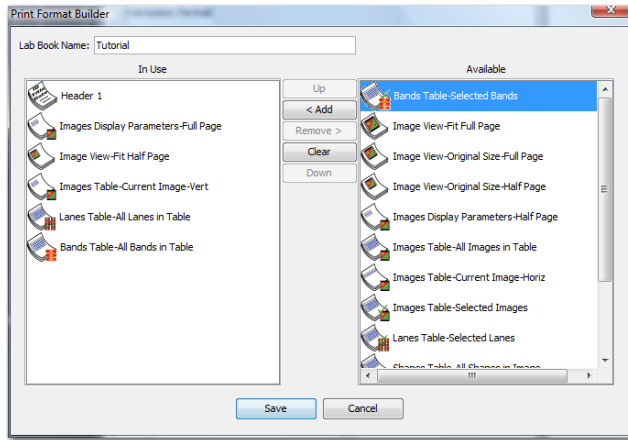

- d. Click once on **Bands Table-Selected Bands** to select it and click **<Add** to add it to the *In Use* column. This Table will now appear in the report.
- e. Click once on **Bands Table-All Bands in Table** to select it and click **Remove>** to take it out of the *In Use* column. This Table will not appear in the report.
- f. Click once on an item in the *In Use* column and click **Up** or **Down** to move the item up or down in the order of the report.
- g. Click **Save** to include the items in the *In Use* column in the report.
- 3) In the **Print** menu, click on **Tutorial** and then **Print** to print the custom report for the current image.